



# User Guide for GoldBio Buffers



**GoldBio © 2015 All Rights Reserved**

**This Publication is a creation of Gold Biotechnology and is intended as a sourcebook for research laboratories. This publication may not be redistributed or copied for commercial use.**

Procedure for use with Gold Biotechnology Buffers;

## Table of Contents

Introduction .....	5
What is a buffer and how do they work? .....	5
Why is this important for biologists? .....	6
<i>Why are there so many different buffers and how do I decide which one to use?</i> .....	6
<i>What are Good's Buffers?</i> .....	6
Table 1. List of Good's Buffers offered by GoldBio in order of pKa. ....	8
How to Choose the Right GoldBio Buffer for Your Experiment .....	9
How to Make Your Buffer Solution .....	10
Temperature Effects .....	11
GoldBio Buffers Grouped By Structural Family .....	12
Acetamido Family (ACES, ADA) .....	12
ACES .....	13
ADA .....	14
Bis(2-hydroxyethyl)amine Family (Bis-Tris, BES, Bicine) .....	15
Bis-Tris .....	15
BES .....	16
Bicine .....	17
Cyclohexylamino Family (CHES, CAPS) .....	17
CHES .....	18
CAPS .....	19
Morpholinic Family (MES, MOPS, MOPSO) .....	19
MES .....	20
MOPS .....	21
MOPSO .....	22
Piperazinic Family (PIPES, HEPES, HEPPSO) .....	22
PIPES .....	23

**Gold Biotechnology**

St. Louis, MO

Ph: (800)248-7609

Web: [www.goldbio.com](http://www.goldbio.com)

Email: [contactgoldbio86@goldbio.com](mailto:contactgoldbio86@goldbio.com)

HEPES .....	24
HEPPSO .....	26
Tris Family (Tris, Bis-Tris, TES, TAPS, Tricine) .....	26
Bis-Tris .....	26
TES .....	27
Tricine .....	27
Tris .....	29
TAPS .....	30
Other buffers .....	30
PBS (Phosphate Buffered Saline) .....	30
Boric Acid .....	31
FAQS .....	32
Troubleshooting .....	33
Protocols .....	34
0.1M ACES Buffer - 1 L .....	35
0.5M ADA Buffer - 1 L .....	36
1M BES Buffer - 1 L .....	37
1M Bicine Buffer - 1 L .....	38
1M Bis-Tris Buffer - 1 L .....	39
0.5M CAPS Buffer - 1 L .....	40
0.5M CHES Buffer - 1 L .....	41
1M HEPES Buffer - 1 L .....	42
1M HEPES-Na Buffer - 1 L .....	43
1M HEPPSO Buffer - 1 L .....	44
0.5M MES Buffer - 1 L .....	45
0.5M MES Monohydrate Buffer - 1 L .....	46
0.5M MES-Na Buffer - 1 L .....	47
1M MOPS Buffer - 1 L .....	48
1M MOPS-Na Buffer - 1 L .....	49
0.5M MOPSO Buffer - 1 L .....	50
0.5M MOPSO-Na Buffer - 1 L .....	51
1M PIPES Buffer - 1 L .....	52
1M PIPES-Na Buffer - 1 L .....	53

1M TAPS Buffer - 1 L .....	54
1M TES Buffer - 1 L.....	55
1M Tricine Buffer - 1 L.....	56
1M Tris Buffer - 1 L.....	57
1M Tris HCl Buffer - 1 L .....	58
Additional Resources .....	59
Biological Buffers Resource Page.....	59
Stock Solution Calculator .....	59
Helpful Articles.....	59
GoldBio Buffers by Category .....	60
Table 2. Biological Buffers.....	60
Table 3. Chromatography Buffers.....	60
Table 4. Cell Culture Buffers .....	61
Table 5. Electrophoresis Buffers .....	61
Table 6. Lysis Buffers.....	61
Biological Buffer Selection Guide.....	62
Biological Buffer Selection Guide (part 2).....	63
Biological Buffer Selection Guide (part 3).....	64
Guide: Commonly used Buffers .....	65
Guide: Commonly used Buffers (part 2) .....	66
Ultimate Buffer Selection Guide .....	67
References .....	68

© 2015 by Gold Biotechnology

All rights reserved. No part of this document may be reproduced or transmitted in any form or by any means, electronic, mechanical, photocopying, recording, or otherwise, without prior written permission of Gold Biotechnology.

Editor – Christopher Harper  
Revised: 10/8/2019

GoldBio · FM-000009 Rev. 2.0  
TD – 1.0  
Creation Date: 9/5/2015


## User Guide for GoldBio Buffers

### Introduction


#### What is a buffer and how do they work?

A buffering agent undergoes readily reversible proton exchange, picking up or dropping off hydrogen ions as they become available. A buffer solution can therefore maintain a stable pH upon the addition of an acid or base more easily than a simple solution of water. Another way to define a buffer is a solution comprised of a weak acid (proton donor, HA) and its conjugate base (proton acceptor, A<sup>-</sup>), or a weak base and its conjugate acid. For instance, a buffer solution can be made of acetic acid and sodium acetate. Such a buffer solution will be between pH 3-6 and will resist changes in pH upon addition of acid or base within that pH range.

HA can dissociate into A<sup>-</sup> and H<sup>+</sup>. H<sup>+</sup> then reacts with water to form H<sub>3</sub>O<sup>+</sup>. In the aqueous buffer solution, H<sub>3</sub>O<sup>+</sup>, HA and H<sup>+</sup> exist in equilibrium with each other. The buffering mechanism consists of two reversible reactions where the concentration of proton donor and proton acceptor are equal.

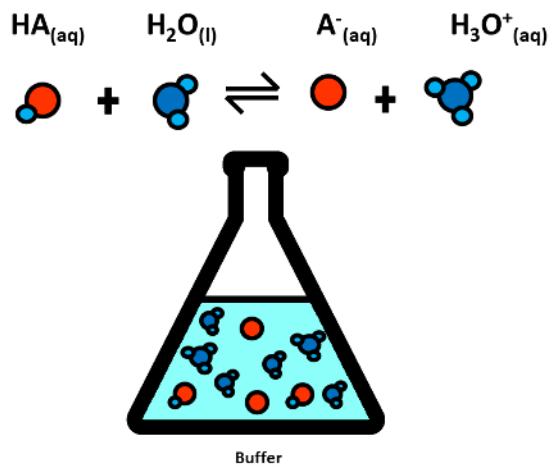

Then, when a strong acid or base is introduced into this system by the scientist or by enzymatic activity during the experiment, the new ions from the introduced acid or base (H<sup>+</sup> or OH<sup>-</sup>) are absorbed by the buffer and the pH remains stable preventing changes in protein structure and function.

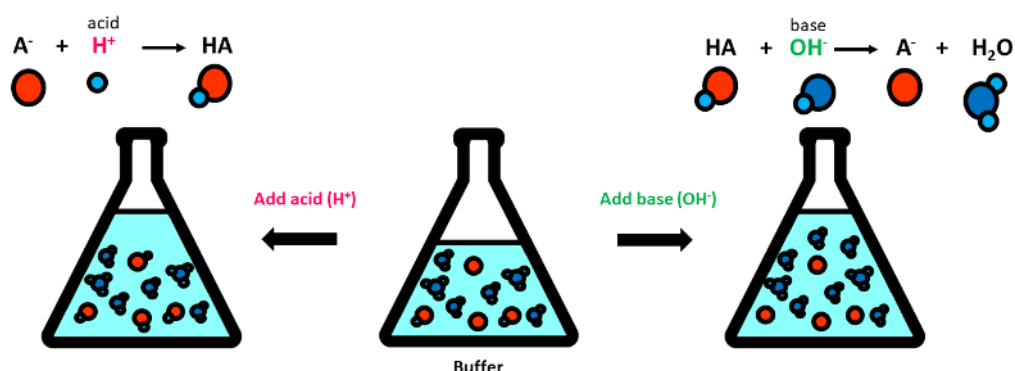

Buffers cannot arbitrarily moderate any changes in ion concentration. Their optimal **buffering capacity, or range**, is defined by the dissociation constant, or  $K_a$ , of the acid. We commonly discuss buffering capacity in terms of the  $pK_a$  or the logarithmic constant of  $K_a$ . We consider the buffering capacity of a specific buffer to be the  $pK_a \pm 1$ . For example, a buffer with a pH of 6.8 has a pH buffering range of 5.8-7.8.

### Why is this important for biologists?

Since protonation alters the overall molecular charge, and can subsequently affect the structure of a molecule, most biological molecules are only functional in a certain pH range, and therefore, many organisms can only survive in a certain pH range. Researchers use buffers in the laboratory to maintain pH conditions that promote stable and functional molecules or organisms.

### Why are there so many different buffers and how do I decide which one to use?

Buffers are chosen based upon the experiment that will be performed. For example, an Acetate buffer will work very well for maintaining a pH between 3 and 6, but it is of little use to you if you're studying infectious agents that spread through blood since blood usually maintains a pH around 7.4. While a phosphate buffer will provide the proper pH range, a phosphate buffer might not be appropriate if the experiment involves an enzyme that is affected by phosphate, such as many kinases, phosphatases, dehydrogenases, carboxypeptidases, ureases and more. Oftentimes, a zwitterionic buffer, such as one of Good's buffers, will work best. There are many different buffers, each with unique, physical properties that allow each individual buffer to have different conditions which are appropriate for use. Sometimes trying to decide which buffer to use can be difficult. Gold Biotechnology, Inc. has created this handbook to aid in the process of choosing a buffer so that you can spend *less* time looking and *more* time doing.

### What are Good's Buffers?

Good's buffers are a set of biochemical buffers chosen by Norman Good based on a set of criteria essential to biological research. In total, there are 10 features that a buffer should

meet in order for it to be useful in research.

1. **pK<sub>a</sub> is between 6 - 8:** Good's criteria were developed in order to identify buffers ideal for biochemical experiments. Therefore, the first criterion is that a biological buffer should be within a neutral pH range, matching biological systems – a pK<sub>a</sub> ranging between 6 and 8.
2. **Soluble in water:** Because biochemical reactions typically take place in aqueous environments, a buffer needs to be water-soluble.
3. **Membrane impermeability:** A biological buffer should not readily cross the cell membrane. Zwitterionic buffers such as MOPS and HEPES do not pass through the cell membrane.
4. **Minimal salt effects:** Since biological buffers are used in biological systems, high salt content could have disastrous effects on your sample. Therefore, buffers should have negligible amounts of ionic content.
5. **Influences on dissociation:** Temperature, concentration and ionic strength can all have an impact on buffers in an experiment. However, for biochemical work, an ideal buffer would be one whose concentration, temperature and ionic composition of the medium will only have minimal effect on buffering capability (pK<sub>a</sub>).
6. **Well-behaved cation interactions:** The formation of complexes between a metal ion and the buffer results in proton release, which affects the pH of the system and may have an adverse effect on experimental results. Thus, these ionic complexes should be soluble and their binding constant must be known. A buffer with a low metal-binding constant is suitable for the study of metal-dependent enzymatic reactions. If your experimental design requires the use of a metal, then you should choose a buffer that does not form a complex with that specific metal.
7. **Stability:** Buffers should be stable and resist enzymatic and nonenzymatic degradation. And they should not interfere with enzyme substrates or resemble them.
8. **Biochemical inertness:** An ideal biological buffer will be inert, meaning that it won't influence or have an active role in biochemical reactions.
9. **Optical absorbance:** Because of the way biological buffers are used, they should not absorb light in the visible or ultraviolet regions of the spectrum. This criterion prevents interference in spectrophometric assays.
10. **Ease of Preparation and Affordability:** It is also important that a biological buffer be easy to use, from preparation to purification. And, they must be affordable.



“[...] IT MAY BE THAT THE QUEST FOR  
UNIVERSAL BIOLOGICAL INERTNESS IS  
FUTILE.”

Norman E. Good

From these criteria, Good *et al.* developed 20 buffers that are commonly referred to as “Good’s buffers.” Many other buffers have been developed to fit these criteria and are often referred to as “Good’s buffers,” despite not being developed by Good himself. No single buffer seems to fulfill all of these criteria, so selecting a buffer for one’s own experiment often involves trial and error before selecting a suitable buffer. In the final analysis, the buffer must fulfill as many criteria as possible and sufficient controls should be designed to exclude unanticipated buffer effects.

The first property of a buffer that should be considered is the *pH range* required for the experiment. Every buffer has a pH range where the buffering capacity is highest, which is generally determined by the  $pK_a$  of the acid (or  $pK_a$ s for polyprotic acids). **Choose a buffer that has  $pK_a$  that is at least within one pH unit of the target pH.** If studying a protein, the target pH will likely be a pH at which the protein is stable, which is generally at least one pH unit away from the *pI* of the protein.

Next, potential deleterious or unwanted buffer *interactions* with native molecules must be considered. Analyze the components of your system and determine, if possible, if there are any potential reactive species. Properly designed controls are critical to rule out buffer effects.

**Table 1. List of Good’s Buffers offered by GoldBio in order of  $pK_a$ .**

Buffer	$pK_a$ at 20°C	$pK_a$ /°C	Catalog Numbers
MES	6.15	-0.011	M-090, M-095, M-091
ADA	6.6	-0.011	A-780
PIPES	6.8	-0.0085	P-280, P-281
ACES	6.9	-0.020	A-010
MOPSO	6.95	-0.015	M-795, M-790
MOPS	7.15	-0.013	M-791, M-790
BES	7.15	-0.016	B-780
TES	7.5	-0.020	T-785
HEPES	7.55	-0.014	H-400, H-401
Tricine	8.15	-0.021	T-870
Tris	8.2		T-400
Bicine	8.35	-0.018	B-785
TAPS	8.55	-0.025	T-780
CHES	9.3		C-870
CAPS	10.4		C-040



## How to Choose the Right GoldBio Buffer for Your Experiment

### Questions to consider:

- *Are there metals in the solution?* Many buffers form a complex with certain metals. If this is the case with your preferred buffer, the concentration of the buffer may need to be adjusted and a stability constant for the buffer-metal complex should be taken into account. You might also want to consider the addition of a metal chelator, such as [EDTA](#).
- *Are there carbohydrates, nucleotides or proteins in the solution?* Borate is known to form a complex with a wide variety of biological molecules. Most amine based buffers that have a neutral pH range, interact with DNA, which can have an effect on electrophoretic mobility and restriction enzyme kinetics.
- *Is your target protein an enzyme and if so, to what class of enzymes does it belong?* Phosphate and Tris, both widely used as buffers, can act as inhibitors for many different enzymes. Other buffers may have unintended, previously unknown effects as well.
- *What are the ionic strength requirements for ligand binding, protein stability or electrophoresis?* The ionic strength of the solution can affect the stability of some enzymes and the reaction kinetics with an enzyme substrate. The ionic strength of the solution will also affect electrophoretic mobility and conductivity during electrophoresis. This is important to keep in mind when preparing a buffer solution. A solution of “free acid” buffer that is brought to its pH with NaOH will have a different ionic strength than the sodium salt of that same buffer brought to the final pH with HCl due to the formation of NaCl in the latter example. The best way to avoid this is to combine the free acid buffer with the sodium salt of that buffer in molar proportions based on the Hendersen-Hasselbalch equation (see [How to Make Your Buffer Solution](#)). Mixing two buffers to achieve the desired pH allows the investigator to avoid the use of strong acids or bases.
- *Are there enzyme inhibitors in the solution?* Enzyme inhibitors are sometimes subject to interactions with buffers. [DEPC](#), a popular RNase inhibitor, is inactivated by Tris and HEPES. However, DEPC is unaffected by MOPS or PBS. If Tris buffer must be used, treat the water with DEPC first, autoclave to destroy the DEPC and then Tris buffer may be used.
- *Does the study involve reduction-oxidation reactions?* HEPES and PIPES have demonstrated the ability to form radical species. Tricine in the presence of peroxide and enzymes with oxidase activity forms a stable nitric oxide-tricine radical species. MES, MOPS, HEPES and Tris have demonstrated the ability to slow Fe<sup>2+</sup> autoxidation.
- *What sort of downstream assays and methods will be used?* Most zwitterionic buffers are known to interfere with the Lowry method of protein quantitation. Tris has been

#### Gold Biotechnology

St. Louis, MO

Ph: (800)248-7609

Web: [www.goldbio.com](http://www.goldbio.com)

Email: [contactgoldbio86@goldbio.com](mailto:contactgoldbio86@goldbio.com)

observed to interfere with the Bradford assay. Some buffers have also been observed to interfere with protein quantitation using Bicinchoninic Acid (BCA). This interference with color development during the BCA assay can sometimes be overcome by precipitating the protein with deoxycholate and trichloroacetic acid or acetone. ADA absorbs UV light at 260 nm which can interfere with spectrophotometric nucleic acid quantitation.

- *Is there any literature regarding interactions of the buffer with any component of the system?* An effective literature search can sometimes go a long way in ruling out disruptive buffers for your experiment.
- Combining two buffers can provide advantages when compared to a single buffer system. While HEPES is a great buffer at physiological pH, the lowering of pH due to metabolic acidification in a cell culture can eventually bring the media out of the desired pH range and cause inhibition of cell growth. Bicarbonate/ CO<sub>2</sub> buffers alone are usually avoided because CO<sub>2</sub> can leave or enter the system freely, unintentionally changing the pH. Waldman *et al.* has shown that creating a buffering system that uses HEPES and bicarbonate can be more effective at maintaining a neutral extracellular pH of a cell culture than either buffer alone. Bicarbonate helps neutralize acid produced during metabolism while HEPES helps maintain the neutral pH after the addition of bicarbonate and after fluctuations of dissolved CO<sub>2</sub> levels.

The last thing to consider is the *concentration* of buffer you need to effectively maintain the pH. If the concentration of buffer is too low, it will not be effective. If the concentration is too high, there exists a greater chance that the buffer will interfere with the experiment. If the system being buffered does not actively exchange hydrogen protons, then a buffer concentration of 25-100mM is usually a good starting point. If proton exchange is likely to occur, then a buffer concentration that is 20X more than the molarity of the protons being exchanged should be used (**NOTE:** this is not always possible to calculate.) The buffer concentration also affects the ionic strength of the solution and hence, the electrophoretic mobility and conductivity during electrophoresis.

## How to Make Your Buffer Solution

For instructions to make a specific buffer solution, see the protocols section at the end of this handbook.

A buffer solution is all about pH!

pH is a way to express the concentration of hydrogen protons in solution. This is relevant because the concentration of hydrogen protons will determine the ratio of protonated to deprotonated species of a biological molecule.

$$\text{pH} = -\log[\text{H}^+]$$

$\text{pK}_a$  is the ionization constant; it is the pH at which a molecule becomes protonated or deprotonated, resulting in equal concentrations of the two chemical species. The buffering capacity is greatest, and therefore, a buffer works best to resist changes in pH, when the pH of a solution is at the  $\text{pK}_a$  value of a buffer. Each buffer should be used in a solution that is within one pH unit of the  $\text{pK}_a$  of the buffer. pH and  $\text{pK}_a$  can be related by the *Henderson-Hasselbalch* equation:

$$\text{pH} = \text{pK}_a + \log\left(\frac{[\text{A}^-]}{[\text{HA}]}\right)$$

Wherein, **[A<sup>-</sup>]** represents the concentration of the **conjugate base** and **[HA]** represents the **concentration of the acid**.

The *Henderson-Hasselbalch* equation can be used to determine the concentration of buffer acid and buffer salt needed to achieve the desired pH. For example, if you are trying to make a 1M solution of HEPES at pH 7.2 (25°C) you can insert the  $\text{pK}_a$  and the desired pH into the Henderson-Hasselbalch equation to find the ratio of HEPES free acid and HEPES sodium salt needed.

$7.20 = 7.48 + \log\left(\frac{[\text{HEPES sodium salt}]}{[\text{HEPES free acid}]}\right)$  This can also be expressed as follows:

$$\frac{[\text{HEPES sodium salt}]}{[\text{HEPES free acid}]} = 0.72$$

This ratio can be divided by two to determine the amount of sodium salt to use so that 0.36mol/L of HEPES sodium salt should be combined with 0.64mol/L of HEPES free acid to obtain a 1M solution with pH 7.2.

Zwitterionic free acids can also be made into a buffer solution using NaOH. Zwitterionic sodium salts can be made into a buffer solution using HCl, but this will result in the formation of NaCl as well.

## Temperature Effects

It is important to know that the  $\text{pK}_a$  of a molecule and the pH of a solution changes when the temperature changes.  $d(\text{pK}_a)/dT$  is an estimate of the change in  $\text{pK}_a$  with respect to temperature. The greater the absolute value of the  $d(\text{pK}_a)/dT$ , the more the  $\text{pK}_a$  will change as the temperature changes. As the  $\text{pK}_a$  changes, the pH of the solution will change and the

buffering capacity will be greatest at the new  $pK_a$ . For some buffers the change is negligible and for others the change is great enough to consider using a different buffer if there will be dramatic temperature changes. For example, when using PIPES or ADA the pH drops 0.24 pH units when the temperature rises from 4°C to 37°C, but the  $pK_a$  and pH of a solution of Tris can change almost an entire pH unit as the temperature changes from 4°C to 37°C.

**When preparing a buffer solution, it is important to make the solution at the temperature and concentration at which you intend to use it.** Dissociation is dependent on temperature and concentration and can sometimes shift dramatically when one of these variables changes, affecting the pH of a solution. If diluting a stock solution of a buffer with high ionic strength, such as PIPES or ADA, be sure to check the pH of your working solution and adjust it if necessary. If the temperature of the media is going to fluctuate widely, be sure to choose a buffer than can accommodate these changes.

**Buffers with the greatest temperature dependence:**

Tris	-0.028
Tricine	-0.021
TES	-0.020
ACES	-0.020
TAPS	+0.018
Bicine	-0.018

## GoldBio Buffers Grouped By Structural Family


Most of the molecules deemed suitable as buffers for biological work can be grouped into six different structural families: Acetamido (ACES, ADA), Bis(2-hydroxyethyl)amine (BES, Bicine, Bis-Tris), Cyclohexylamino (CAPS, CHES), Morpholinic (MES, MOPS, MOPSO), Piperazinic (HEPES, HEPPSO, PIPES), and Tris (Tris, TES, TAPS, Tricine, Bis-Tris).

### Acetamido Family (ACES, ADA)

The Acetamido family includes [ACES](#) and [ADA](#). Both of these buffering agents are known to complex with various metals. Despite the stronger acidity of the sulfonic acid group on ACES, the presence of the less substituted amine gives ACES a slightly higher  $pK_a$  than the diacetic amine substitution imparts to ADA.

## ACES

pH range: 6.1 – 7.5

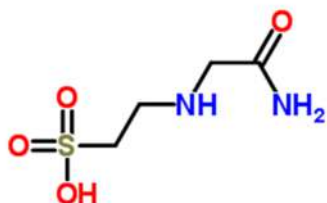

GoldBio Catalog: [A-010](#)

IUPAC Name: 2-[[2-amino-2-oxoethyl]amino]ethanesulfonic acid

Formula: C<sub>4</sub>H<sub>10</sub>N<sub>2</sub>O<sub>4</sub>S

MW: 182.20 g/mol

pK<sub>a</sub> at 25°C: 6.78

Solubility at 0°C: 0.22M


**Metal complexation:** strong complex with Mg<sup>2+</sup> and Cu<sup>2+</sup>;  
weak complex with other common metals

ACES is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. ACES is capable of forming complexes with some common metals, so formation constants should be taken into account when using this buffer in a solution containing metal ions. ACES absorbs UV light at 230 nm, which may interfere with some spectrophotometric assays. ACES is used as buffering component in cell culture media, in both capillary and gel electrophoresis (Liu *et al.* 1999, Hamoudová *et al.* 2006), and in various protein studies. It has been used as an enzyme assay buffer for the study of glucosidases (Hoskins *et al.* 1985), as a buffer for x-ray crystallography of aldehyde dehydrogenase enzyme complexes (Perez-Miller *et al.* 2003), and as a buffer solution to wash and heat *Lactobacillus plantarum* cells for the purpose of studying the effects of heat stress on the duration of the lag phase of growth of such cells (Smelt *et al.* 2002). It is used in the preparation of buffered charcoal yeast extract agar for the isolation of *Legionella* spp. (Edelstein *et al.* 1993), as well as in buffered culture media for hairy roots of *Catharanthus roseus* (Morgan *et al.* 2000). It has been observed to act as a competitive inhibitor of GABA receptor binding to rat brain synaptic membranes **and its use as a buffer should thus be avoided when studying GABA receptor interactions** (Tunnicliff *et al.* 1981).

**ACES Keywords:** *Legionella*, BCYE agar, x-ray crystallography, electrophoresis, *Catharanthus roseus*, GABA inhibition, glucosidases

## ADA

**pH range: 6.0 – 7.2**

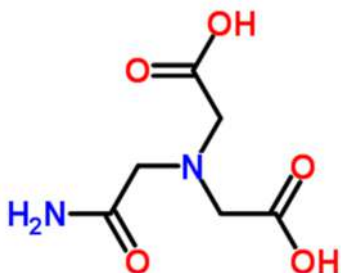

**GoldBio Catalog:** [A-780](#)

**IUPAC Name:** 2,2'-[(2-Amino-2-oxoethyl)imino]diacetic acid

**Formula:** C<sub>6</sub>H<sub>10</sub>N<sub>2</sub>O<sub>5</sub>

**MW:** 190.22 g/mol

**pK<sub>a</sub> at 25°C:** 6.59

**Solubility at 0°C:** solubility of free acid is very small; solubility of sodium salts very large

**Metal complexation:** strong complex with Mn<sup>2+</sup>, Co<sup>2+</sup>, Ni<sup>2+</sup>, Zn<sup>2+</sup>, Cd<sup>2+</sup>, Pb<sup>2+</sup> and Cu<sup>2+</sup>; weak complex with other common metals

ADA is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. ADA is practically insoluble in water, but solubility of ADA increases if the free acid is converted to the sodium salt with NaOH (see [ADA Stock Solution protocol](#)). **ADA forms a strong complex with many common metals so stability constants should be taken into account when using this buffer.** While ADA can be used as a buffer in cation exchange chromatography, it should be used in lower concentrations because of the relatively large ionic strength and dependence of concentration on pK<sub>a</sub>. ADA absorbs UV light between 0.1 and 260 nm, which may interfere with spectrophotometric assays.

**ADA will not dissolve until you add a strong base!**


ADA buffer has been used successfully as a solution for protein crystallization (Wilson 2003, Dahl *et al.* 2007) and as a solution that permits microtubule polymerization to occur (Weisenberg *et al.* 1972). It was used as a running buffer during capillary electrophoresis and determined to provide the best resolution of the five buffers tested when separating amiodarone and desethylamiodarone (Pérez-Ruiz *et al.* 2002). ADA has been used as a sample buffer during differential scanning calorimetry to study [FGF1](#), which could not be studied in phosphate or sulfate based buffers (Blaber *et al.* 1999). ADA has also been used as a complexing agent to effectively help remove contaminant metals from soil (Chen *et al.* 1995). ADA is not suitable for use with the Bicinchoninic Acid (BCA) Assay.

**ADA Keywords:** crystallization, running buffer, capillary electrophoresis, FGF-1, differential scanning calorimetry

## Bis(2-hydroxyethyl)amine Family (Bis-Tris, BES, Bicine)

The Bis(2-hydroxyethyl)amine family includes Bis-Tris, BES, and Bicine. These buffering agents are known to complex with various metals. The substituted groups on the main moiety are, in order of increasing acidity: a carboxylic acid group ([Bicine](#)), a sulfonic acid group ([BES](#)), and a Tris(hydroxymethyl) group ([Bis-Tris](#)).

### Bis-Tris

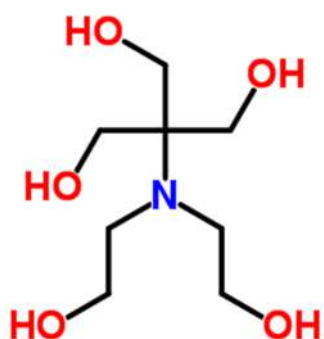

pH range: 5.8 – 7.2

GoldBio Catalog: [B-020](#)

IUPAC Name: 2-[Bis(2-hydroxyethyl)amino]-2-(hydroxymethyl)-1,3-propanediol

Formula: C<sub>8</sub>H<sub>19</sub>NO<sub>5</sub>

MW: 209.24 g/mol

pK<sub>a</sub> at 25°C: 6.46

Solubility at 0°C: >1M

**Metal complexation:** strong complex with Pb<sup>2+</sup> and Cu<sup>2+</sup>;

weak complex with other common metals

Bis-Tris, also known as Bis-Tris Methane, is a zwitterionic buffering agent that contains functional groups related to two different structural families, the Bis(2-hydroxyethyl)amine family and the Tris family. Bis-Tris forms a strong complex with Cu<sup>2+</sup> and Pb<sup>2+</sup> and a weak complex with many common metals, so formation constants should be taken into account when using this buffer in a solution containing metal ions. Bis-Tris has been suggested as a substitute for the highly toxic buffer cacodylate (Scopes 1994).


Try replacing cacodylate with a safer alternative like bis-tris for crystallization.

Bis-Tris buffered solutions are useful for resisting changes in pH during pressure changes (Funtenberger *et al.* 1995). Bis-Tris is used as a sample buffer, gel buffer and running buffer with various types of electrophoresis (Nijtmans *et al.* 2002, Klampfl *et al.* 1998) and as a buffer during anion exchange chromatography (Donoghue *et al.* 2000). It has also been used during x-ray crystallography of chloroplastic oxygen evolving enhancer protein (Kohoutova *et al.* 2009). Bis-Tris has been determined to be a useful buffer for NMR spectroscopy because of the low conductivity and high sensitivity that can be achieved using it (Kelly *et al.* 2002). Bis-Tris has been reported to interact with human liver fatty acid binding protein (FAB) and affect protein dynamics (Long *et al.* 2009). Bis-Tris is not suitable for use with the Bicinchoninic Acid (BCA) Assay.

**Bis-Tris Keywords:** electrophoresis, anion exchange chromatography, NMR spectroscopy, cacodylate

## BES

**pH range: 6.4 – 7.8**

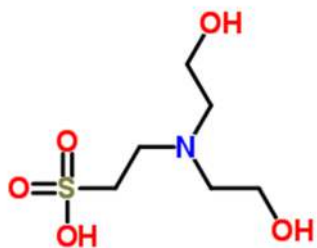

**GoldBio Catalog:** [B-780](#)

**IUPAC Name:** 2-[Bis(2-hydroxyethyl)amino]ethanesulfonic acid

**Formula:** C<sub>6</sub>H<sub>15</sub>NO<sub>5</sub>S

**MW:** 213.25 g/mol

**pK<sub>a</sub> at 25°C:** 7.09

**Solubility at 0°C:** 3.2M

**Metal complexation:** weak complex with Co<sup>2+</sup> and Cu<sup>2+</sup>


BES is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. BES can be used as a binding buffer and eluent in cation exchange chromatography as well as a buffer in gel filtration chromatography (Schirch *et al.* 1985). BES has been used in culture media as a buffer during bacteriophage adsorption (Lu *et al.* 2003). In an attempt to culture embryonic chondrocytes, it was found that using BES resulted in the adverse formation of cytoplasmic vacuoles, which were absent in the non-zwitterionic buffered media (Poole *et al.* 1982). BES forms a complex with DNA, affecting restriction enzyme kinetics. It interferes with reactions between DNA and restriction enzymes to a lesser extent than similar buffers with less substituted amine groups, such as Tris, but to a greater extent than HEPES (Wenner *et al.* 1999). BES can be used in a buffered saline solution in the calcium phosphate mediated transfection of eukaryotic cells with plasmid DNA (Chen *et al.* 1987, Schenborn 2000). Though BES binds to copper ions, it is still suitable for use with the Bicinchoninic Acid (BCA) Assay.

**BES, like most neutral buffers that contain an amine group, forms a complex with DNA that affects restriction enzyme kinetics.**

**BES Keywords:** cation exchange chromatography, gel filtration, bacteriophage adsorption, calcium phosphate transfection, BCA assay



## Bicine

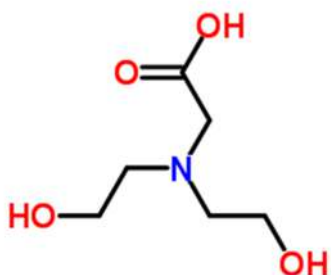

**pH range: 7.6 – 9.0**

**GoldBio Catalog:** [B-785](#)

**IUPAC Name:** N,N-Bis(2-hydroxyethyl)glycine

**Formula:** C<sub>6</sub>H<sub>13</sub>NO<sub>4</sub>

**MW:** 163.17 g/mol

**pK<sub>a</sub> at 25°C:** 8.26

**Solubility at 0°C:** 1.1M

**Metal complexation:** strong complex with Mg<sup>2+</sup>, Ca<sup>2+</sup>, Fe<sup>3+</sup>, Co<sup>2+</sup> and Cu<sup>2+</sup>; weak complex with Mn<sup>2+</sup>

Bicine is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. Bicine forms a strong complex with many common metals; therefore, stability constants and concentrations should be taken into consideration when choosing this buffer. It has been used in solutions for protein crystallization (Kanyo *et al.* 1992) and as a mobile phase buffer and eluent in cation exchange chromatography. Bicine can be used as a running buffer in doubled SDS-PAGE and has been reported to give better resolution and representation of proteins than the use of glycine or tricine (Williams *et al.* 2006). Bicine is commonly used as a buffer in the polymerase chain reaction, serving as both a pH buffer and a metal ion buffer (Innis *et al.* 1999). It has also been used in a multiphasic buffer system for SDS-PAGE of proteins (Wiltfang *et al.* 1991). Bicine has been used as an extraction/grinding buffer for extraction of enzymes from various plant species, including salt-sensitive plants (Greenway *et al.* 1972, Yeoh *et al.* 1981). It was determined to be a suitable buffer for the culture of ammonia fungi at pH 8 (Yamanaka 2003). Bicine has been determined to be a useful buffer for NMR spectroscopy because of the low conductivity and high sensitivity that can be achieved using it (Kelly *et al.* 2002). Bicine is not suitable for use with the Bicinchoninic Acid (BCA) Assay.

**For the best resolution during doubled SDS-PAGE, try using bicine.**


**Bicine Keywords:** crystallization, cation exchange chromatography, salt-sensitive plants, ammonia fungi, NMR spectroscopy, dSDS-PAGE, PCR

## **Cyclohexylamino Family (CHES, CAPS)**

The Cyclohexylamino family includes [CHES](#) and [CAPS](#). **Neither buffer is known to form any significant complex with metals.** They are used for studies occurring above physiological pH. The longer carbon chain of the sulfonic acid group in CAPS gives it a slightly more alkaline range than CHES.

## CHES

**pH range: 8.6 – 10.0**

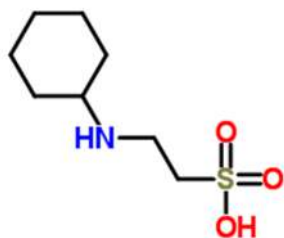

**GoldBio Catalog:** [C-870](#)

**IUPAC Name:** 2-(Cyclohexylamino)ethanesulfonic acid

**Formula:** C<sub>8</sub>H<sub>17</sub>NO<sub>3</sub>S

**MW:** 207.29 g/mol

**pK<sub>a</sub> at 25°C:** 9.5

**Solubility at 0°C:** 1.14M


**Metal complexation:** no significant complexes observed

CHES is a zwitterionic buffering agent used in biochemistry and molecular biology that is sometimes referred to as a Good's buffer despite not being introduced by Good *et al.* It is commonly used to study enzymatic processes above physiological pH. CHES displays only weak or negligible complexation capabilities for some metals making it a suitable noncoordinating buffer for use in most solutions that contain metal ions. It has been used as a buffer for enzymatic reactions (Omburo *et al.* 1992, Wasserman *et al.* 1984) and as a sample and running buffer for various types of capillary electrophoresis (Marina *et al.* 1996, Lavigne *et al.* 2007, Cunliffe *et al.* 2002). CHES displays high affinity for the iodoacetate binding site of liver alcohol dehydrogenase, protecting the enzyme from inactivation by iodoacetic acid (Syvertsen *et al.* 1981). Teresa Thiel *et al.* (1998) observed the greatest phosphodiesterase activity in CHES compared with the other buffers tested and showed that CHES can inhibit the viability of *Micrococcus luteus* at pH 9.5.

**If your experiment requires alkaline conditions, consider using CHES or CAPS.**

**CHES Keywords:** alkaline, capillary electrophoresis, phosphodiesterase, noncoordinating

## CAPS

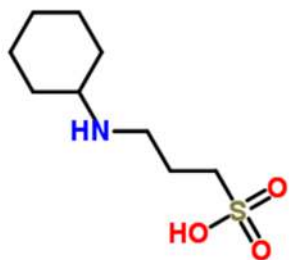

**pH range: 9.7 – 11.1**

**GoldBio Catalog:** [C-040](#)

**IUPAC Name:** 3-(Cyclohexylamino)-1-propanesulfonic acid

**Formula:** C<sub>9</sub>H<sub>19</sub>NO<sub>3</sub>S

**MW:** 221.32 g/mol

**pK<sub>a</sub> at 25°C:** 10.4

**Solubility at 0°C:** 0.47M

**Metal complexation:** no significant complexes observed


CAPS is another zwitterionic buffering agent used in biochemistry and molecular biology that was not introduced by Good *et al.*, but is sometimes referred to as a Good's buffer. It is commonly used to study enzymatic processes above physiological pH. CAPS displays only weak or negligible complexation capabilities for some metals, making it a suitable noncoordinating buffer for use in solutions that contain metal ions. CAPS has been used as a transfer buffer for diffusion blotting and electroblotting (Thomas *et al.* 1996, Antonetti *et al.* 1999, Hulett *et al.* 1990), as a binding buffer and eluent in cation exchange chromatography (Thomas *et al.* 1996), as a running buffer in capillary electrophoresis (Prange *et al.* 1999), as an effective crystallization solution for various proteins (Zaitseva *et al.* 2004, Celie *et al.* 2005), and as an enzyme assay buffer (Hulett *et al.* 1990, Donarski *et al.* 1989, Jennings *et al.* 2001). CAPS can support alkaline phosphatase activity and inhibit the growth of *Aeromonas* spp. ATCC 31085 at pH 10.5 (Thiel *et al.* 1998). CAPS is suitable for use with the Bicinchoninic Acid (BCA) Assay.

**CAPS Keywords:** transfer buffer, Western blot, electroblotting, cation exchange, alkaline phosphatase, BCA

## **Morpholinic Family (MES, MOPS, MOPSO)**

The Morpholinic family includes [MES](#), [MOPS](#), and [MOPSO](#). These buffers are generally ideal for use in solutions that contain metal ions. The longer carbon chain of the sulfonic acid in MOPS makes it slightly less acidic than MES, while the additional hydroxyl group in MOPSO makes it slightly more acidic than MOPS.

## MES

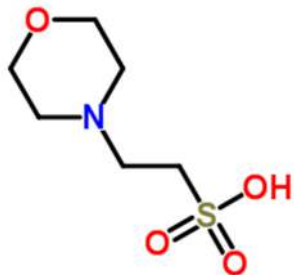

**pH range: 5.5 – 6.7**

**GoldBio Catalog:** [M-090](#), [M-095](#), [M-091](#)

**IUPAC Name:** 2-(4-Morpholinyl)ethanesulfonic acid

**Formula:** C<sub>6</sub>H<sub>13</sub>NO<sub>4</sub>S

**MW:** 195.24 g/mol

**pK<sub>a</sub> at 25°C:** 6.1

**Solubility at 0°C:** 0.65M

**Metal complexation:** possible complex with Fe<sup>3+</sup>

MES is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. MES lacks the ability to form a complex with most metal ions and is recommended for use as a noncoordinating buffer in solutions with metal ions. MES is sometimes employed as a substitute for the highly toxic cacodylate, as well as a substitute for the ionic buffers citrate and malate. MES is often used in buffered culture media for bacteria, yeast and mammalian cells. It is toxic towards most plants at high concentrations, but can be used in plant culture media at concentrations of ~10mM (Parfitt *et al.* 1988).


Looking for a nontoxic, zwitterionic alternative to cacodylate, citrate, or malate buffers? Try using MES.

Due to low ionic mobility and low conductivity at high concentrations, MES is regarded as an excellent buffer for use in capillary electrochromatography (Boughtflower *et al.* 1995). Those same properties make it **useful in many types of electrophoreses and chromatography** including gel-filtration chromatography (Weisenberg *et al.* 1995), phosphocellulose column chromatography (Alonso *et al.* 2000), hydrophobic interaction chromatography (Alonso *et al.* 2000), cation exchange chromatography (Cook *et al.* 1994), and SDS-PAGE (Kashino *et al.* 2001). MES has seen use been used for fluorescence microscopy (Neill *et al.* 2002), as an assay buffer to study Tau protein (Alonso *et al.* 2000), in a culture medium to initiate growth of pine and fir trees (Pullman *et al.* 2005), and in culture media for the growth of halophilic bacteria (Rodriguez-Valera *et al.* 1980). Heteromeric Cx32/Cx26 connexin channels isolated from mouse liver were seen to be inhibited by MES at any pH near or below its pK<sub>a</sub> (Bevans *et al.* 1999). MES has been observed to complex with some organotin(IV) molecules that are being explored as antitumor agents (Mohamed *et al.* 2007) and to bind human liver fatty acid binding protein which affected the protein dynamics (Long *et al.* 2009).

**MES Keywords:** culture media, chromatography, running buffer, electrophoresis, fluorescence microscopy, Tau

## MOPS

**pH range: 6.5 – 7.9**

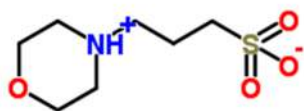

**GoldBio Catalog:** [M-790](#), [M-791](#)

**IUPAC Name:** 3-(Morpholin-4-ium-4-yl)-1-propanesulfonate

**Formula:** C<sub>7</sub>H<sub>15</sub>NO<sub>4</sub>S

**MW:** 209.26 g/mol

**pK<sub>a</sub> at 25°C:** 7.14

**Solubility at 0°C:** large

**Metal complexation:** possible complex with Fe<sup>3+</sup>

MOPS is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. MOPS lacks the ability to form a complex with most metal ions and is recommended for use as a noncoordinating buffer in solutions with metal ions. MOPS is often used in cell culture media for bacteria, yeast and mammalian cells; though, concentrations higher than 20mM are not recommended for mammalian cell work. It is commonly used for cell culture media, as a running buffer in electrophoresis and for protein purification in chromatography.


MOPS has been employed in cell culture media for the growth of enterobacteria (Neidhardt *et al.* 1974) and *Lactococcus lactis* (Jensen *et al.* 1993), in culture media to study meiotic regulation in mouse oocytes (Downs *et al.* 1997), and as an incubation buffer for MAP kinases isolated from rat cells (Zhang *et al.* 1997). In a study characterizing an atrazine metabolizing enzyme, de Souza *et al.* (1996) used it as a lysis buffer for *E. coli* cells, as an enzyme assay buffer for atrazine chlorohydrolase, as an eluent in gel filtration chromatography, and as a buffer during mass spectrometry. MOPS has been used to measure absorption during UV/VIS spectrophotometry and to study redox properties using cyclic voltammetry (Collins *et al.* 2005). MOPS is regarded as an excellent buffer for use in separating RNA in agarose gels, due to the high resolutions separation that occurs and a lack of interaction between MOPS and the RNase-inhibitor DEPC (Brody *et al.* 2004). It has also seen use as a running buffer in SDS-PAGE (Reichel *et al.* 2009). It is recommended to sterilize MOPS buffers by filtration rather than with autoclave due to the unknown identity of yellow degradation products that occur after sterilization of MOPS with autoclave. It is suitable for use in the Bicinchoninic Acid (BCA) Assay.

**Do not autoclave MOPS!  
Sterilize by filtration to avoid  
degradation.**

**MOPS Keywords:** culture media, running buffer, electrophoresis, chromatography, RNA, SDS-PAGE, BCA assay

### MOPSO

**pH range: 6.2 – 7.6**

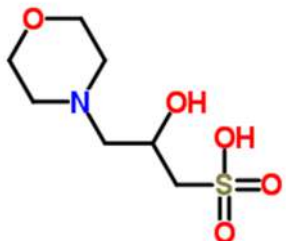

**GoldBio Catalog:** [M-795](#), [M-799](#)

**IUPAC Name:** 2-Hydroxy-3-(4-morpholinyl)-1-propanesulfonic acid

**Formula:** C<sub>7</sub>H<sub>15</sub>NO<sub>5</sub>S

**MW:** 225.26 g/mol

**pK<sub>a</sub> at 25°C:** 6.87

**Solubility at 0°C:** 0.75M

**Metal complexation:** possible complex with Fe<sup>3+</sup>


MOPSO is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. MOPSO lacks the ability to form a complex with most metal ions and is recommended for use as a noncoordinating buffer in solutions with metal ions. MOPSO has been used as a carrier electrolyte in capillary electrophoresis (Zemann *et al.* 2001), as a grinding buffer for the extraction of sesquiterpene synthases from *Artemisia annua* (Bouwmester *et al.* 1999), and as a crystallization buffer for homoglutathione synthetase (Galant *et al.* 2009). It has been used during fluorescence spectroscopy, spectrophotometry, and isothermal titration calorimetry (Ukaegbu *et al.* 2010). MOPSO has also been successfully used to maintain the pH in solutions of raw sugar, thereby preventing color fluctuations (Ahmedna *et al.* 1997). Colby *et al.* used it as a lysis buffer and enzyme assay buffer in a study characterizing spearmint limonene synthase from recombinant *E. coli* (1993). MOPSO interacts with the peptide backbone of [bovine serum albumin \(BSA\)](#) to stabilize BSA against thermal denaturation (Taha *et al.* 2011).

**MOPSO Keywords:** capillary electrophoresis, crystallization, fluorescence spectroscopy, isothermal titration Calorimetry (ITC), bovine serum albumin

### **Piperazinic Family (PIPES, HEPES, HEPPSO)**

The Piperazinic family includes [PIPES](#), [HEPES](#) and [HEPPSO](#). With the exception of a HEPPSO-Cu<sup>2+</sup> complex, these buffers do not generally form complexes with metals. **They are generally not suitable for reduction-oxidation studies.** The ethanol group in place of one of the ethanesulfonic acid groups makes HEPES more alkaline than PIPES. The electron donating hydroxyl group present on the ethanesulfonic group of HEPPSO gives it a more alkaline nature than HEPES.

## PIPES

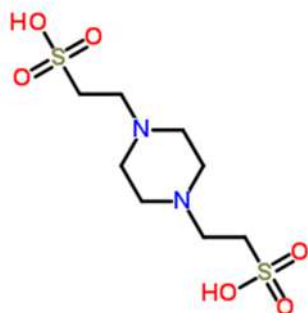

**pH range: 6.1 – 7.5**

**GoldBio Catalog:** [P-281](#), [P-280](#)

**IUPAC Name:** 2,2'-(1,4-Piperazinediyl)diethanesulfonic acid

**Formula:** C<sub>8</sub>H<sub>18</sub>N<sub>2</sub>O<sub>6</sub>S<sub>2</sub>

**MW:** 302.37 g/mol

**pK<sub>a</sub> at 25°C:** 6.76

**Solubility at 0°C:** solubility of free acid is very small; solubility of sodium salts very large

**Metal complexation:** no significant complexes observed

PIPES is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. **PIPES is practically insoluble in water, but solubility of PIPES increases if the free acid is converted to the sodium salt with NaOH** (see [PIPES Stock Solution protocol](#)). PIPES lacks the ability to form a significant complex with most metal ions and is recommended for use as a noncoordinating buffer in solutions with metal ions. This buffer is capable of forming radicals and is therefore not suitable for redox reactions (Baker *et al.* 2007). While **PIPES can be used as a buffer in cation exchange chromatography, it should be used in lower concentrations** because of the relatively large ionic strength and dependence of concentration on pK<sub>a</sub>.


PIPES is regarded as an excellent buffer for the fixation of plant cells for electron microscopy (Bauer *et al.* 1977), and a suitable replacement for the highly toxic cacodylate buffer in this application (Yamamoto *et al.* 1983). It has been used to purify microtubule proteins using phosphocellulose chromatography (Clayton *et al.* 1980), in [PHEM](#) buffer alongside HEPES to study cytoskeletal components (Schliwa *et al.* 1981), to purify recombinant GTP-binding proteins, ARF1 and ARF2, by gel filtration (Cockcroft *et al.* 1994), and as a buffer solution to crystallize a transketolase enzyme from *E. coli* (Littlechild *et al.* 1995). PIPES has been observed to interact with a truncated variant of human bile salt-stimulate lipase (Moore *et al.* 2001) as well as inhibit a K<sup>+</sup>-independent phosphatase reaction catalyzed by a (Na<sup>+</sup> + K<sup>+</sup>) –ATPase from dog kidney (Robinson *et al.* 1987). It is suitable for use in the Bicinchoninic Acid (BCA) Assay.

**PIPES is an excellent buffer to use in place of cacodylate buffer for electron microscopy.**

**PIPES Keywords:** radicals, electron microscopy, plant cell fixation, microtubules, PHEM, crystallization, BCA assay, cell culture



## HEPES

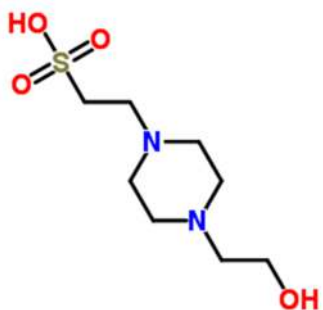

**pH range: 6.8 – 8.2**

**GoldBio Catalog:** [H-400](#), [H-401](#)

**IUPAC Name:** 2-[4-(2-Hydroxyethyl)-1-piperazinyl]ethanesulfonic acid

**Formula:** C<sub>8</sub>H<sub>18</sub>N<sub>2</sub>O<sub>4</sub>S

**MW:** 238.30 g/mol

**pK<sub>a</sub> at 25°C:** 7.48

**Solubility at 0°C:** 2.25M

**Metal complexation:** no significant complexes observed

HEPES is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. It meets many of Good's criteria for an effective buffer. It is useful at physiological pH, it does not readily complex with any metals, it is very soluble in water, it is relatively cheap, and it is biologically inert in most cases. An exception to its inertness is the participation in reduction-oxidation reactions to generate free radicals (Grady *et al.* 1988, Kirsch *et al.* 1998, Yang *et al.* 1999, Baker *et al.* 2007). HEPES is commonly used in cell culture media for various types of organisms and for protein studies that use it as a binding buffer and an eluent during cation exchange chromatography (Yan *et al.* 1998, Susin *et al.* 1999). It has been used to study pH effects on seed germination (Basto *et al.* 2013) and as a grinding buffer in plant studies (Soltis *et al.* 1983), as a running buffer (with or without Tris) in various forms of gel electrophoresis (Michels *et al.* 2002, Koort *et al.* 2002, Praveen 2009), and to flatten the pH gradient in polyacrylamide gels containing pH 5-7 ampholines during ultrathin-layer isoelectric focusing (Budowle 1987). It is not uncommon for HEPES to be used in DNA studies such as a buffer solution to form precipitates of calcium phosphate and DNA (Wigler *et al.* 1979), as a deposition buffer for atomic force microscopy (Lyubchenko *et al.* 2011), and as a buffer in electroporation (Fromm *et al.* 1985). HEPES interferes with reactions between DNA and restriction enzymes, but to a lesser extent than similar buffers with less substituted amine groups, such as Tris (Wenner *et al.* 1999). This lack of interference occurs mostly because steric hindrance of the protonated tertiary amine in HEPES allows for a weaker association with the negative charge of DNA than most

HEPES, like most neutral buffers that contain an amine group, forms a complex with DNA that affects restriction enzyme kinetics. Due to steric hindrance, this effect is minimal with HEPES when compared with Tris.



buffers, shielding DNA from restriction enzymes to a lesser extent than less substituted buffers.

Itagaki and Kimura (1979) demonstrated that HEPES can be used to buffer mammalian cell cultures in closed systems of high and low density cultures, as well as open systems of high density cultures. Zigler *et al.* (1985) showed that HEPES was capable of increasing the phototoxicity of mammalian tissue culture media containing riboflavin in a study done with murine thymocytes. This phototoxicity is attributed to the inhibition of thymidine uptake caused by the generation of hydrogen peroxide. Photoreactivity may be decreased by keeping HEPES-buffered media in the dark or by lowering or eliminating the amount of riboflavin, tryptophan and tyrosine present in media (Stoien *et al.* 1974). A fair amount of studies have demonstrated the usefulness of HEPES as buffer for *in vitro* fertilization and embryo culture (Mahadevan *et al.* 1986, Behr *et al.* 1990)). **Heteromeric Cx32/Cx26 connexin channels isolated from mouse liver were seen to be inhibited by HEPES at any pH near or below its pK<sub>a</sub>, indicating that HEPES is a poor buffer to use for studying such heteromeric connexin channels** (Bevans *et al.* 1999). The research of Habib and Tabata (2004) led to an understanding of the free radicals generated by HEPES when Au<sup>3+</sup> was present. Their research indicated that the damage to DNA was caused by nitrogen-centered cationic free radicals and not reactive oxygen species (ROS). Supporting the addition of HEPES to Tris-based buffers, Ahmad and Ghasemi (2007) showed that **adding HEPES to the traditional Tris buffer used in quantitative real-time PCR provided lower background levels and expanded dynamic range**. HEPES has been deemed suitable for use as a buffer in quantitative protein measurement with the Bicinchoninic Acid assay (BCA) (Kaushal *et al.* 1986), but unsuitable for use with the Lowry (Folin) method of protein quantitation (Peterson 1979).

**HEPES Keywords:** cation exchange, running buffer, electrophoresis, chromatography, DNA, electroporation, phototoxicity with riboflavin, *in vitro* fertilization (IVF), quantitative real time polymerase chain reaction (qPCR), restriction enzymes, atomic force microscopy, seed germination, radicals

## HEPPSO

**pH range: 7.1 – 8.5**

**GoldBio Catalog:** [H-480](#)

**IUPAC Name:** 2-Hydroxy-3-[4-(2-hydroxyethyl)-1-piperazinyl]-1-propanesulfonic acid


**Formula:** C<sub>9</sub>H<sub>20</sub>N<sub>2</sub>O<sub>5</sub>S

**MW:** 268.33 g/mol

**pK<sub>a</sub> at 25°C:** 7.84

**Solubility at 0°C:** 2.20M

**Metal complexation:** weak complex with Cu<sup>2+</sup>

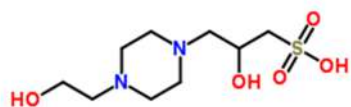

HEPPSO is a zwitterionic buffer that was selected and described by Good *et al.* in 1966. HEPPSO has been used as a crystallization buffer solution for glycogen synthase enzymes and for enzymes complexed with HEPPSO (Sheng *et al.* 2009). HEPPSO displayed little to no toxicity for the indicator species *Daphnia magna* and has thus been deemed suitable for toxicological studies (Keating *et al.* 1996). This buffer is capable of forming radicals and is therefore not suitable for redox reactions. HEPPSO forms a complex with Cu<sup>2+</sup> ions; therefore, stability constants and concentrations should be taken into consideration when using this buffer in a solution containing copper ions. Though HEPPSO binds to copper ions, it is still suitable for use with the Bicinchoninic Acid (BCA) Assay (Kaushal *et al.* 1986).

**HEPPSO is suitable for toxicity studies.**

**HEPPSO Keywords:** crystallization, glycogen synthase, nontoxic, BCA assay, radicals

## **Tris Family (Tris, Bis-Tris, TES, TAPS, Tricine)**


The Tris family includes [Tris](#), [Bis-Tris](#), [TES](#), [TAPS](#), and [tricine](#). These buffers are known to form complexes with various metals. They are widely used in electrophoresis.

### Bis-Tris

See description under [Bis\(2-hydroxyethyl\)amine Family](#) heading.

## TES

**pH range: 6.8 – 8.2**

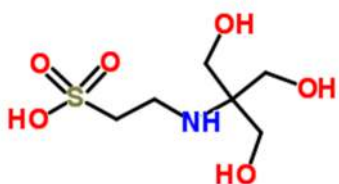

**GoldBio Catalog:** [T-785](#)

**IUPAC Name:** 2-[[1,3-Dihydroxy-2-(hydroxymethyl)-2-propanyl]amino]ethanesulfonic acid

**Formula:** C<sub>6</sub>H<sub>15</sub>NO<sub>6</sub>S

**MW:** 229.25 g/mol

**pK<sub>a</sub> at 25°C:** 7.40

**Solubility at 0°C:** 2.60M


**Metal complexation:** strong complex with Cr<sup>3+</sup> and Fe<sup>3+</sup>; weak complex with Co<sup>2+</sup>, Ni<sup>2+</sup>, Cu<sup>2+</sup>, and Zn<sup>2+</sup>

TES is a zwitterionic buffering agent used in biochemistry and molecular biology that was selected and described by Good *et al.* It has a **pK<sub>a</sub> that is right at physiological pH (7.4) at 25°C**, making it a useful buffer for many biologists. Comai & Kosuge (1980) used TES as a cell wash buffer and as a solution from which to precipitate and extract DNA. TES has also been used to help purify microsomes from porcine kidney by gel filtration and affinity chromatography, and was used in the same study as an enzyme assay buffer (Mumford *et al.* 1981). TES is used in Test Yolk Buffer, which is known as a first choice solution for preserving spermatozoa (Weidel *et al.* 1987, Hammadeh *et al.* 2001). Like most neutral pH amine-based buffers, TES forms a complex with DNA; it affects restriction enzymes kinetics more than HEPES but less than Tris (Wenner *et al.* 1999). TES has been used in media for viral plaque assays (Richter 1967). It is suitable for use in the Bicinchoninic Acid (BCA) Assay.

**TES Keywords:** physiological pH, affinity chromatography, gel filtration, Test Yolk Buffer, viral plaque assay, BCA assay

## Tricine

**pH range: 7.4 – 8.8**

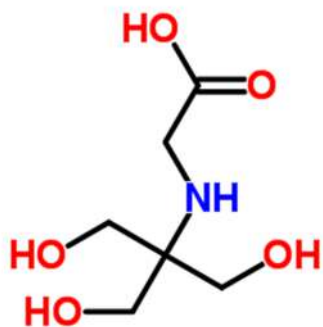

**GoldBio Catalog:** [T-870](#)

**IUPAC Name:** N-[1,3-Dihydroxy-2-(hydroxymethyl)-2-propanyl]glycine

**Formula:** C<sub>6</sub>H<sub>13</sub>NO<sub>5</sub>

**MW:** 179.17 g/mol

**pK<sub>a</sub> at 25°C:** 8.05

**Solubility at 0°C:** 0.8M

**Metal complexation:** strong complex with  $Mg^{2+}$ ,  $Ca^{2+}$ ,  $Ni^{2+}$ ,  $Co^{2+}$ ,  $Zn^{2+}$  and  $Cu^{2+}$

Tricine is a zwitterionic buffering agent used in biochemistry and molecular biology that was selected and described by Good *et al.* Tricine is commonly used instead of [glycine](#) in the Laemmli buffer during SDS-PAGE to separate small peptides (Schägger 1987). It is used in animal tissue culture (Gardner 1969) and helps inhibit the unintended growth of mycoplasma (Spendlove *et al.* 1971). Tricine has also been added to bacterial culture media to prevent precipitation of iron salts (Neidhardt *et al.* 1974). Tricine was determined to allow more light production from firefly luciferase in the presence of ATP than other buffers tested (Webster *et al.* 1980). Tricine has been used as an enzyme assay buffer for isocitrate dehydrogenase and glutamate dehydrogenase (Gambord *et al.* 1970).


Try using tricine in place of the controlled substance barbital when following protocols that use barbital as a buffer.

Tricine can bind hydroxyl radicals, forming formaldehyde in the process (Shiraishi *et al.* 1993, Hicks *et al.* 1986). It can also form tricine nitroxide radicals; this should be taken into consideration if using tricine in the presence of an enzyme with oxidase activity (Grande *et al.* 1978). **Tricine forms a strong complex with many common metals so stability constants and concentrations should be taken into consideration when choosing this buffer.** Even though tricine forms a complex with  $Cu^{2+}$ , the addition of  $Cu^{2+}$  to a solution enables tricine to be used in the Lowry protein assay. Tricine is not suitable for use with the Bicinchoninic Acid (BCA) Assay. Tricine has been used as a substitute for protocols that utilize the controlled substance barbital as a buffer for electrophoresis (Monthony *et al.* 1978).

**Tricine Keywords:** Laemmli buffer, SDS-PAGE, animal tissue culture, bacterial culture media, firefly luciferase, isocitrate dehydrogenase, glutamate dehydrogenase, radicals, Lowry protein assay, barbital

## Tris

pH range: 7.0 – 9.0

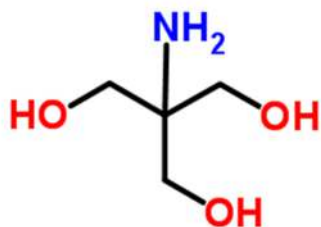

GoldBio Catalog: [T-400](#), [T-095](#), [T-090](#)

IUPAC Name: 2-Amino-2-(hydroxymethyl)-1,3-propanediol

Formula: C<sub>4</sub>H<sub>11</sub>NO<sub>3</sub>

MW: 121.14 g/mol

pK<sub>a</sub> at 25°C: 8.06

Solubility at 0°C: 2.40M


**Metal complexation:** strong complex with Cr<sup>3+</sup>, Fe<sup>3+</sup>, Ni<sup>2+</sup>, Co<sup>2+</sup>, and Cu<sup>2+</sup>; weak complex with other common metals

Short for Tris (hydroxymethyl) aminomethane (THAM), Tris is an organic compound often used in buffer solutions for gel electrophoresis, such as TAE or TBE. It is used in the preparation of Laemmli buffer, one of the most common running buffers for SDS-PAGE. In addition, Tris can also be used for many custom running and loading buffers, most often with [glycine](#) and SDS. It can be used as a buffer and eluent in anion exchange chromatography (Huisman *et al.* 1965). The use of **single-junction pH electrodes that contain silver should be avoided** when determining the pH of a solution containing Tris buffer (Benesch *et al.* 1955). **The pH of a solution containing Tris is very dependent on temperature; while temperature and concentration should be taken into to consideration for every buffer, it is especially important to prepare Tris buffer at the temperature at which it will be used** (El-Harakany *et al.* 1984, Vega *et al.* 1985). The primary amine group of Tris reacts, to some extent, with a variety of molecules, including the RNase inhibitor [DEPC](#), many aldehydes, common metals, many enzymes and DNA. This reactivity combined with the tendency of Tris to permeate membranes **excludes it from most cell culture work**. Tris is not suitable for use with the Bicinchoninic acid (BCA) assay.

**Do not use single-junction pH electrodes in Tris buffer.**

**Tris Keywords:** running buffers, electrophoresis, TAE, TBE, Laemmli, glycine, SDS-PAGE, anion exchange chromatography, temperature dependent

## TAPS

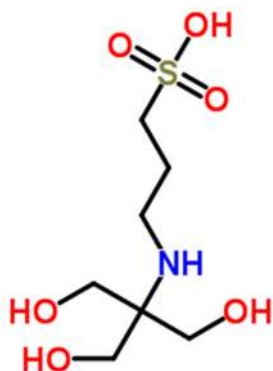

**pH range: 7.7 – 9.1**

**GoldBio Catalog:** [T-780](#)

**IUPAC Name:** 3-[[1,3-Dihydroxy-2-(hydroxymethyl)-2-propanyl]amino]-1-propanesulfonic acid

**Formula:** C<sub>7</sub>H<sub>17</sub>NO<sub>6</sub>S

**MW:** 243.28 g/mol

**pK<sub>a</sub> at 25°C:** 8.40

**Solubility at 0°C:** large

**Metal complexation:** strong complex with Cr<sup>3+</sup>, Fe<sup>3+</sup>, and Cu<sup>2+</sup>; weak complex with other common metals

TAPS is a zwitterionic buffering agent used in biochemistry and molecular biology that was selected and described by Good *et al.* TAPS has been used in planar chromatography to separate dyes (Nurok *et al.* 2000), in ion exchange chromatography to purify glutathione S-transferase isolated from atrazine-resistant *Abutilon theophrasti* (Anderson *et al.* 1991), and as an enzyme assay buffer for aminopeptidase activity (Nishiwaki *et al.* 2000). It is a preferred culture media buffer to use for dinoflagellate experiments, allowing for minimal pH change and maximal growth (Loeblich *et al.* 1975). Heteromeric Cx32/Cx26 connexin channels isolated from mouse liver were seen to be inhibited by TAPS at any pH near or below its pK<sub>a</sub> (Bevans *et al.* 1999). TAPS forms a strong complex with Cr<sup>3+</sup>, Fe<sup>3+</sup>, and Cu<sup>2+</sup> and a weak complex with some other common metals, so stability constants and concentrations should be taken into account when choosing this buffer.

**TAPS Keywords:** planar chromatography, ion exchange chromatography, glutathione S-transferase, dinoflagellate, aminopeptidase, connexin inhibition

## Other buffers

### PBS (Phosphate Buffered Saline)

**pH range: 7.2 – 7.4**

**GoldBio Catalog:** [P-271](#)


PBS (phosphate buffered saline) is a common buffer solution used in the laboratory. [PBS buffer tablets from GoldBio](#) make a solution that contains 10mM phosphate buffer, 137mM sodium chloride, and 2.7mM potassium chloride. It is an isotonic and nontoxic

buffer that is meant to mimic the pH, osmolarity, and ion concentrations of the human body, making it particularly useful in tissue cultures at physiological pH (7.4). PBS is used for immunoassays, for sample dilution, for protein purification, and as a wash buffer in cell cultures. The addition of [EDTA](#) to PBS allows the solution to be used to disengage attached or clumped cells because of the chelation of  $\text{Ca}^{2+}$  and  $\text{Mg}^{2+}$  ions involved in adhesion.

**PBS Keywords:** isotonic, physiological pH, immunoassays, sample dilution, cell culture

### Boric Acid

**pH range: 8.2 – 10.1**

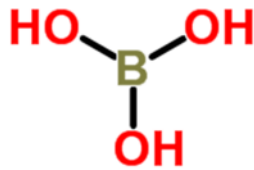

**GoldBio Catalog:** [B-030](#), [B-031](#)

**IUPAC Name:** Boric acid

**Formula:**  $\text{H}_3\text{BO}_3$

**MW:** 61.83 g/mol

**pK<sub>a</sub> at 25°C:** 9.23

**Solubility at 0°C:** 0.4M

Boric acid is commonly used for the preparation of buffers such as [Tris-Borate-EDTA](#) (TBE). Borate buffers have been used in capillary and gel electrophoresis for separating proteins, nucleic acids, and secondary metabolites such as catecholamines (Kaneta *et al* 1991). The useful buffering range of boric acid buffers is pH 8.2 - 10.1. Boric acid and its related salts form a complex equilibria of at least 10 reactions, and at pH 8 or higher, the tetraborate ion ( $\text{B}_4\text{O}_7^{2-}$ ) is the most predominant species present (Salentine, *et al.* 1983, Trejo *et al.* 2012). Boric acid is used in catalytic amounts to promote amidation between carboxylic acids and amines (Tang 2005). It is capable of forming a complex with various carbohydrates (Aronoff *et al.* 1975), as well as with pyridine nucleotides, such as [NAD<sup>+</sup>](#) (Johnson *et al.* 1976). Boron is also an essential micronutrient for plants and can be supplied in the form of boric acid (Warington 1923).

**Boric Acid Keywords:** TBE, electrophoresis, plant micronutrient, amidation

## FAQS

### ***Why are the solubilities given at 0°C if I'm going to be making my buffer at room temperature?***

Norman Good determined the solubility of all the buffers with which he worked at 0°C. This is important to know if the buffer will be used in an ice bath or cold room so that precipitation of the buffer can be avoided.

### ***Why won't ADA or PIPES go into solution?***

Both ADA and PIPES have very limited solubility in their free acid forms. Once conditions become less acidic, usually through the addition of NaOH, these buffers will go into solution.

### ***It doesn't look like Bicine or Tricine is going into solution. What do I do?***

Add the full volume of water and mix well to make [Bicine](#) or [Tricine](#) at the concentration suggested in each protocol. **NOTE:** A lower concentration should be used if the solution is to be frozen.

### ***What's the difference between Tris-HCl and Tris base?***

Tris HCl is used to simplify the process of making Tris buffer solutions. It also allows for more reproducibility. Rather than adding HCl or NaOH to adjust the pH, you can use Tris-HCl instead. You can find [more details in this article](#).

### ***My buffer solutions look a little cloudy. Is this okay?***

Is the buffer in question MOPS? MOPS will sometimes have a light yellow appearance. Swirl the bottle to see if anything settles at the bottom or if the solution remains consistent. If it is dark yellow or if the solution is not consistent, dispose of the solution properly. For other buffers, a cloudy appearance is likely occurring due to contamination of microbes. It is also possible that your buffer solution has become chemically unstable. Dispose of the solution properly and prepare new buffer solutions.

### ***How do I make an RNase-free solution containing Tris or HEPES?***

The reactive amine group on Tris reacts with diethylpyrocarbonate (DEPC); therefore, when doing RNA work be sure you do not treat Tris-containing solutions with DEPC. HEPES also interferes with the inactivation of RNases by DEPC. You can treat a solution of deionized water with DEPC to inactivate any RNases, autoclave the solution to inactivate the DEPC, and then add the buffer. Instead, consider using [GoldBio MOPS](#) for RNA work.

### ***How do I choose a buffer in which my protein will be soluble and active?***



Make sure that the pH for which you're aiming is far away from the isoelectric point (pI) of your protein. Then, read to find out what other researchers before you have used on similar proteins. Try experimenting with different buffers that are in the proper pH range. Ionic strength is very important so try changing the concentration of salt in your buffer solution. Determine if there is a metal needed that is critical for protein structure. Determine if there is any metal in solution that might disrupt the protein structure and if so, try adding a metal chelator like [EDTA](#).

## Troubleshooting

### *Why are my cells dying in HEPES-buffered media?*

If your media contains riboflavin, then HEPES is most likely increasing the phototoxicity of the media to your cells. Try lowering the amount of riboflavin, tyrosine or tryptophan in the media, or try protecting the culture from light by wrapping it in foil.

### *I've used a certain buffer to grow these same embryos before and seen great expression of gene A but now I'm seeing poor expression of gene B. Could my choice of buffer be affecting this observation?*

Gene expression of embryos or oocytes grown *in vitro* can actually be affected by buffer choice (Palasz *et al* 2008).

## Protocols

Protocols to make a buffer solution using each of the GoldBio buffering agents are given in alphabetical order.

### Stock Solutions:

[ACES](#)

[ADA](#)

[BES](#)

[Bicine](#)

[Bis-Tris](#)

[CAPS](#)

[CHES](#)

[HEPES](#)

[HEPES, Sodium](#)

[HEPPSO](#)

[MES](#)

[MES, Monohydrate](#)

[MES, Sodium](#)

[MOPS](#)

[MOPS, Sodium](#)

[MOPSO](#)

[MOPSO, Sodium](#)

[PIPES](#)

[PIPES, Sodium](#)

[TAPS](#)

[TES](#)

[Tricine](#)

[Tris](#)

[Tris HCl](#)

## 0.1M ACES Buffer - 1 L

### Instructions

1. Dissolve 18.22 g of ACES ([ACES, GoldBio Catalog # A-010](#) [CAS 7365-82-4, mw. = 182.20 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 0.1M ACES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 4.09  
Adjust pH with: 10N NaOH

pH	6.1	6.2	6.3	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5
mL	1.1	1.2	1.42	1.72	2.02	2.5	2.9	3.3	3.75	4.25	4.71	5.11	5.49	5.89	6.21

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

ACES pKa at 25°C: 6.78  
ACES pH range: 6.1 -7.2  
d(pKa)/dT value: -0.02

## 0.5M ADA Buffer - 1 L

### Instructions

1. Suspend 95.11 g of ADA ([ADA, GoldBio Catalog # A-780](#) [CAS 26239-55-4, mw. = 190.22]) in 750 mL of dH<sub>2</sub>O. ADA will not readily dissolve until the pH is raised.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 0.5M ADA, use the table below to estimate the required volume of base for a given pH:

Starting pH: 1.95  
Adjust pH with: 10N NaOH

pH	6.0	6.1	6.2	6.3	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2
mL	6.3	6.5	6.75	7	7.25	7.5	7.8	8.1	8.35	8.5	8.7	8.9	9.05

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

ADA pKa at 25°C: 6.59  
ADA pH range: 6.0 – 7.2  
d(pKa)/dT value: -0.011

## 1M BES Buffer - 1 L

### Instructions

1. Dissolve 213.25 g of BES ([BES, GoldBio Catalog # B-780](#) [CAS 10191-18-1, mw. = 213.25 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M BES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 2.55  
Adjust pH with: 10N NaOH

pH	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8
mL	11	12.5	15	19	22	27	32	38	43	50	56.5	63	71	76	82

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

BES pKa at 25°C: 7.09  
BES pH range: 6.4 – 7.8  
d(pKa)/dT value: -0.016

## 1M Bicine Buffer - 1 L

### Instructions

1. Dissolve 163.17 g of Bicine ([Bicine, GoldBio Catalog # B-785](#) [CAS 150-25-4, mw. = 163.17 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M Bicine, use the table below to estimate the required volume of base for a given pH:

Starting pH: 5.30  
Adjust pH with: 10N NaOH

pH	7.6	7.7	7.8	7.9	8.0	8.1	8.2	8.3	8.4	8.5	8.6	8.7	8.8	8.9	9.0
mL	9	11	12	17	21	26	31	36	40	46	52	58	63	69	74

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

Bicine pKa at 25°C: 8.26  
Bicine pH range: 7.6 – 9.0  
d(pKa)/dT value: -0.018

## 1M Bis-Tris Buffer - 1 L

### Instructions

1. Dissolve 209.24 g of Bis-Tris ([Bis-Tris, GoldBio Catalog # B-020](#) [CAS 6976-37-0, mw. = 209.24 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M Bis-Tris, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 9.90  
Adjust pH with: conc. HCl

pH	5.5	5.6	5.7	5.8	5.9	6.0	6.1	6.2	6.3	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2
mL	77	76	74	72	70	67	64	61	57	53	49	44	39	35	30	26	22	18

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

Bis-Tris pKa at 25°C: 6.50  
Bis-Tris pH range: 5.8 – 7.2  
d(pKa)/dT value: -0.017

## 0.5M CAPS Buffer - 1 L

### Instructions

1. Dissolve 110.66 g of CAPS ([CAPS, GoldBio Catalog # C-040](#) [CAS 1135-40-6, mw. = 221.32 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 0.5M CAPS, use the table below to estimate the required volume of base for a given pH:

Starting pH: 5.03  
Adjust pH with: 10N NaOH

pH	9.7	9.8	9.9	10.0	10.1	10.2	10.3	10.4	10.5	10.6	10.7	10.8	10.9	11.0	11.1
mL	6	8	9	11	13	16	18	21	235	26	29	32	34	37	39

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

CAPS pKa at 25°C: 10.4  
CAPS pH range: 9.7 – 11.1  
d(pKa)/dT value: -0.009



## 0.5M CHES Buffer - 1 L

### Instructions

1. Dissolve 103.65 g of CHES ([CHES, GoldBio Catalog # C-870](#) [CAS 103-47-9, mw. = 207.29 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 0.5M CHES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 4.13  
Adjust pH with: 10N NaOH

pH	8.6	8.7	8.8	8.9	9.0	9.1	9.2	9.3	9.4	9.5	9.6	9.7	9.8	9.9	10.0
mL	5	7	11	15	18	20	22	24	27	30	33	35	37	39	41

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

CHES pKa at 25°C: 9.5  
CHES pH range: 8.6 – 10.0  
d(pKa)/dT value: -0.011

## 1M HEPES Buffer - 1 L

### Instructions

1. Dissolve 238.30 g of HEPES ([HEPES, Free Acid, GoldBio Catalog # H-400](#) [CAS 7365-45-9, mw. = 238.30]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of HEPES free acid and HEPES sodium salt ([HEPES, Sodium Salt, GoldBio Catalog # H-401](#) [CAS 75277-39-3, mw. = 260.29]) can be mixed to attain a pH of ~7.5. The pH can be adjusted by increasing the molar ratio of HEPES free acid (more acidic) or HEPES sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M HEPES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 5.63  
Adjust pH with: 10N NaOH

pH	6.8	6.9	7	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8	8.1	8.2
mL	6	8	10	14	17	20	24	28	33	38	44	49	55	61	67

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

HEPES pKa at 25°C: 7.50  
HEPES pH range: 6.8 – 8.2  
d(pKa)/dT value: -0.014

## 1M HEPES-Na Buffer - 1 L

### Instructions

1. Dissolve 260.29 g of HEPES, Sodium Salt ([HEPES, Sodium Salt, GoldBio Catalog # H-401](#) [CAS 75277-39-3, mw. = 260.29]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl. Note: This method will produce about 0.25M – 0.9M NaCl in the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

**Note:** Alternatively, equimolar concentrations of HEPES free acid ([HEPES, Free Acid, GoldBio Catalog # H-400](#) [CAS 7365-45-9, mw. = 238.30]) and HEPES sodium salt can be mixed to attain a pH of ~7.5. The pH can be adjusted by increasing the molar ratio of HEPES free acid (more acidic) or HEPES sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M HEPES, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 10.56  
Adjust pH with: conc. HCl

pH	6.8	6.9	7	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8	8.1	8.2	8.3
mL	75	73	71	68	65	62	59	55	50	46	43	37	32	28	24	21

**Note:** This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

HEPES pKa at 25°C: 7.50  
HEPES pH range: 6.8 – 8.2  
d(pKa)/dT value: -0.014

## 1M HEPPSO Buffer - 1 L

### Instructions

1. Dissolve 268.33 g of HEPPSO ([HEPPSO, GoldBio Catalog # H-480](#) [CAS 68399-78-0, mw. = 268.33]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M HEPPSO, use the table below to estimate the required volume of base for a given pH:

Starting pH: 5.89  
Adjust pH with: 10N NaOH

pH	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8	8.1	8.2	8.3	8.4	8.5
mL	3.9	5.0	6.4	8.2	10.7	13.3	16.2	19.5	24.3	28.3	33.5	38.8	43.3	49.0	56.0

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

HEPPSO pKa at 25°C: 7.85  
HEPPSO pH range: 7.1 – 8.5  
d(pKa)/dT value: -0.010

## 0.5M MES Buffer - 1 L

### Instructions

1. Suspend 97.62 g of MES free acid ([MES, Free Acid, GoldBio Catalog # M-095](#) [CAS 4432-31-9, mw. = 195.24 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

**Note:** Alternatively, equimolar concentrations of MES free acid and MES sodium salt ([MES, Sodium Salt, GoldBio Catalog # M-091](#) [CAS 71119-23-8, mw. = 217.22]) can be mixed to attain a pH of ~ 6.1. The pH can be adjusted by increasing the molar ratio of MES free acid (more acidic) or MES sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 0.5M MES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 3.23  
Adjust pH with: 10N NaOH

pH	5.5	5.6	5.7	5.8	5.9	6	6.1	6.2	6.3	6.4	6.5	6.6	6.7
mL	5	6.2	7.6	9.3	11.4	13.6	15.8	18.3	21.1	23.7	26.4	28.7	31.1

**Note:** This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MES pKa at 25°C: 6.10  
MES pH range: 5.5 – 6.7  
d(pKa)/dT value: -0.011

## 0.5M MES Monohydrate Buffer - 1 L

### Instructions

1. Suspend 106.53 g of MES free acid monohydrate ([MES, Free Acid Monohydrate, GoldBio Catalog # M-090](#) [CAS 145224-94-8, mw. = 213.25 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

**Note:** Alternatively, equimolar concentrations of MES free acid monohydrate and MES sodium salt ([MES, Sodium Salt, GoldBio Catalog # M-091](#) [CAS 71119-23-8, mw. = 217.22]) can be mixed to attain a pH of ~ 6.1. The pH can be adjusted by increasing the molar ratio of MES free acid (more acidic) or MES sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 0.5M MES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 2.92  
Adjust pH with: 10N NaOH

pH	5.5	5.6	5.7	5.8	5.9	6	6.1	6.2	6.3	6.4	6.5	6.6	6.7
mL	5	6.1	7.4	8.7	10.8	12.6	14.6	16.8	19.2	21.7	24	26.3	28.2

**Note:** This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MES pKa at 25°C: 6.10  
MES pH range: 5.5 – 6.7  
d(pKa)/dT value: -0.011

## 0.5M MES-Na Buffer - 1 L

### Instructions

1. Suspend 108.61 g of MES sodium salt ([MES, Sodium Salt, GoldBio Catalog # M-091](#) [CAS 71119-23-8, mw. = 217.22]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl. This method will produce 0.12M -0.38M NaCl in the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of MES sodium salt and MES free acid ([MES, Free Acid, GoldBio Catalog # M-095](#) [CAS 4432-31-9, mw. = 195.24 g/mol]) can be mixed to attain a pH of ~ 6.1. The pH can be adjusted by increasing the molar ratio of MES free acid (more acidic) or MES sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 0.5M MES, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 10.08  
Adjust pH with: conc. HCl

pH	5.5	5.6	5.7	5.8	5.9	6.0	6.1	6.2	6.3	6.4	6.5	6.6	6.7	6.8
mL	32	31	30	29	28	26	24	22	20	18	16	14	12	10

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MES pKa at 25°C: 6.10  
MES pH range: 5.5 – 6.7  
d(pKa)/dT value: -0.011

## 1M MOPS Buffer - 1 L

### Instructions

1. Dissolve 209.26 g of MOPS free acid ([MOPS Free Acid, Ultra Pure, GoldBio Catalog # M-790](#) [CAS 1132-61-2, mw. = 209.26 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of MOPS free acid and MOPS sodium salt ([MOPS, Sodium Salt, GoldBio Catalog # M-791](#) [CAS 71119-22-7, mw. = 231.25]) can be mixed to attain a pH of ~ 7.2. The pH can be adjusted by increasing the molar ratio of MOPS free acid (more acidic) or MOPS sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M MOPS, use the table below to estimate the required volume of base for a given pH:

Starting pH: 3.82  
Adjust pH with: 10N NaOH

pH	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9
mL	7	9	11	13	16	20	24	28	33	39	45	51	57	62	68

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MOPS pKa at 25°C: 7.14  
MOPS pH range: 6.5 – 7.9  
d(pKa)/dT value: -0.011



## 1M MOPS-Na Buffer - 1 L

### Instructions

1. Dissolve 231.25 g of MOPS, Sodium Salt ([MOPS, Sodium Salt, GoldBio Catalog # M-791](#) [CAS 71119-22-7, mw. = 231.25]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl. Note: This method will produce about 0.24M – 0.88M NaCl in the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of MOPS sodium salt and MOPS free acid ([MOPS Free Acid, Ultra Pure, GoldBio Catalog # M-790](#) [CAS 1132-61-2, mw. = 209.26 g/mol]) can be mixed to attain a pH of ~ 7.2. The pH can be adjusted by increasing the molar ratio of MOPS free acid (more acidic) or MOPS sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M MOPS, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 10.32  
Adjust pH with: conc. HCl

pH	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9
mL	73	70.5	68	66	63	59	55	51	47	42	37	32	28	24	20

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MOPS pKa at 25°C: 7.14  
MOPS pH range: 6.5 – 7.9  
d(pKa)/dT value: -0.011

## 0.5M MOPSO Buffer - 1 L

### Instructions

1. Dissolve 112.63 g of MOPSO, Free Acid ([MOPSO, Free Acid, GoldBio Catalog # M-795](#) [CAS 68399-77-9, mw. = 225.26]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

**Note:** Alternatively, equimolar concentrations of MOPSO free acid and MOPSO sodium salt ([MOPSO, Sodium Salt, GoldBio Catalog # M-799](#) [CAS 79803-73-9, mw. = 247.24]) can be mixed to attain a pH of ~ 6.9. The pH can be adjusted by increasing the molar ratio of MOPSO free acid (more acidic) or MOPSO sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 0.5M MOPSO, use the table below to estimate the required volume of base for a given pH:

Starting pH: 3.47  
Adjust pH with: 10N NaOH

pH	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9
mL	8	10	12	14	16	19	22	25	28	31	33	35	37	39	41

**Note:** This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MOPSO pKa at 25°C: 6.87  
MOPSO pH range: 6.2 – 7.6  
d(pKa)/dT value: -0.015

## 0.5M MOPSO-Na Buffer - 1 L

### Instructions

1. Dissolve 123.62 g of MOPSO sodium salt ([MOPSO, Sodium Salt, GoldBio Catalog # M-799](#) [CAS 79803-73-9, mw. = 247.24]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl. Note: This method will produce about 0.1M – 0.46M NaCl in the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of MOPSO free acid ([MOPSO, Free Acid, GoldBio Catalog # M-795](#) [CAS 68399-77-9, mw. = 225.26]) and MOPSO sodium salt can be mixed to attain a pH of ~ 6.9. The pH can be adjusted by increasing the molar ratio of MOPSO free acid (more acidic) or MOPSO sodium salt (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 0.5M MOPSO, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 10.35  
Adjust pH with: conc. HCl

pH	6.2	6.3	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6
mL	30.8	29.5	28.5	27.3	26	24.4	22.7	20.8	19	17.1	15.1	13.2	11.5	9.6	8.3

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

MOPSO pKa at 25°C: 6.87  
MOPSO pH range: 6.2 – 7.6  
d(pKa)/dT value: -0.015

## 1M PIPES Buffer - 1 L

### Instructions

1. Add 302.37 g of PIPES ([PIPES, Free Acid, GoldBio Catalog # P-281](#) [CAS 5625-37-6, mw. = 302.37]) to 600 mL of dH<sub>2</sub>O. PIPES free acid will not readily dissolve until the solution is raised above pH 6.5.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M PIPES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 2.87  
Adjust pH with: 10N NaOH

pH	6.1	6.2	6.3	6.4	*6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5
mL	78	93	108	110	112	115	119	123	127	132	138	144	148	153	158

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

PIPES pKa at 25°C: 6.76  
PIPES pH range: 6.1 – 7.5  
d(pKa)/dT value: -0.0085

## 1M PIPES-Na Buffer - 1 L

### Instructions

1. Add 335.34 g of PIPES, Sodium Salt ([PIPES, Sodium Salt, GoldBio Catalog # P-280](#) [CAS 100037-69-2, mw. = 335.34]) to 750 mL of dH<sub>2</sub>O. The solution should be at pH ~7.2.
2. Adjust to desired pH using 10N NaOH or concentrated HCl. Using HCl to lower the pH will produce 0.044M – 0.44M NaCl in the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M PIPES, use the table below to estimate the required volume of acid or base for a given pH:

Starting pH: 7.2  
Adjust pH with: 10N NaOH or conc. HCl

<b>pH</b>	6.1	6.2	6.3	6.4	6.5	6.6	6.7	6.8	6.9	7.0	7.1	7.3	7.4	7.5
<b>mL</b>	37.2	35.5	34.0	31.4	29.3	25.4	21.6	17.5	13.3	8.2	3.7	10	15	20
	<u>Concentrated HCl</u>											<u>10N NaOH</u>		
<p><b>Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.</b></p>														

PIPES pKa at 25°C: 6.76  
PIPES pH range: 6.1 – 7.5  
d(pKa)/dT value: -0.0085

## 1M TAPS Buffer - 1 L

### Instructions

1. Dissolve 243.28 g of TAPS ([TAPS, GoldBio Catalog # T-780](#) [CAS 29915-38-6, mw. = 243.28 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M TAPS, use the table below to estimate the required volume of base for a given pH:

Starting pH: 2.81  
Adjust pH with: 10N NaOH

pH	7.7	7.8	7.9	8.0	8.1	8.2	8.3	8.4	8.5	8.6	8.7	8.8	8.9	9.0	9.1
mL	6.5	8	10	13	16	19	23	26	32	36	41	47	52	57	62

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

TAPS pKa at 25°C: 8.40  
TAPS pH range: 7.7 – 9.1  
d(pKa)/dT value: 0.018

## 1M TES Buffer - 1 L

### Instructions

1. Dissolve 229.25 g of TES ([TES, GoldBio Catalog # T-785](#) [CAS 7365-44-8, mw. = 229.25 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M TES, use the table below to estimate the required volume of base for a given pH:

Starting pH: 3.83  
Adjust pH with: 10N NaOH

pH	6.8	6.9	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8.0	8.1	8.2
mL	8	10	13	15	18	22	27	31	36	41	47	52	57	62	67

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

TES pKa at 25°C: 7.40  
TES pH range: 6.8 – 8.2  
d(pKa)/dT value: -0.02

## 1M Tricine Buffer - 1 L

### Instructions

1. Dissolve 179.17 g of Tricine ([Tricine, GoldBio Catalog # T-870](#) [CAS 5704-04-1, mw. = 179.17 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

To make a 1 L solution of 1M Tricine, use the table below to estimate the required volume of base for a given pH:

Starting pH: 5.38  
Adjust pH with: 10N NaOH

pH	7.4	7.5	7.6	7.7	7.8	7.9	8.0	8.1	8.2	8.3	8.4	8.5	8.6	8.7	8.8
mL	9	11	14	17	20	24	28	33	38	43	48	54	59	64	69

**Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.**

Tricine pKa at 25°C: 8.05  
Tricine pH range: 7.4 – 8.8  
d(pKa)/dT value: -0.021



## 1M Tris Buffer - 1 L

### Instructions

1. Dissolve 121.14 g of Tris (Tris Base) ([Tris, GoldBio Catalog # T-400](#) [CAS 77-86-1, mw. = 121.14 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using concentrated HCl.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

**Note:** Alternatively, equimolar concentrations of Tris base and Tris HCl ([Tris HCl, GoldBio Catalog # T-095](#) [CAS 1185-53-1, mw. = 157.60]) can be mixed to attain a pH of ~ 8.1. The pH can be adjusted by increasing the molar ratio of Tris HCl (more acidic) or Tris base (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M Tris, use the table below to estimate the required volume of acid for a given pH:

Starting pH: 10.82  
Adjust pH with: conc. HCl

pH	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8.0	8.1	8.2	8.3	8.4	8.5	8.6	8.7	8.8	8.9	9.0
mL	77	75	73	72	70	67	64	61	57	53	49	45	41	37	32	28	25	21	18	15	12

**Note:** This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

Tris pKa at 25°C: 8.06  
Tris pH range: 7.0 – 9.0  
d(pKa)/dT value: -0.028

## 1M Tris HCl Buffer - 1 L

### Instructions

1. Dissolve 157.60 g of Tris HCl ([Tris HCl, GoldBio Catalog # T-095](#) [CAS 1185-53-1, mw. = 157.60 g/mol]) in 750 mL of dH<sub>2</sub>O.
2. Adjust to desired pH using 10N NaOH. This method will introduce 0.06M - 0.81M NaCl into the concentrated stock solution.
3. Fill to final volume of 1 L with dH<sub>2</sub>O.
4. Filter sterilize (recommended) or autoclave.
5. Store at 4°C.

Note: Alternatively, equimolar concentrations of Tris base ([Tris, GoldBio Catalog # T-400](#) [CAS 77-86-1, mw. = 121.14 g/mol]) and Tris HCl can be mixed to attain a pH of ~ 8.1. The pH can be adjusted by increasing the molar ratio of Tris HCl (more acidic) or Tris base (more basic) and estimated using the Hendersen-Hasselbalch equation.

To make a 1 L solution of 1M Tris HCl, use the table below to estimate the required volume of base for a given pH:

Starting pH: 3.58  
Adjust pH with: 10N NaOH

pH	7.0	7.1	7.2	7.3	7.4	7.5	7.6	7.7	7.8	7.9	8.0	8.1	8.2	8.3	8.4	8.5	8.6	8.7	8.8	8.9	9.0
mL	6	7	9	11	13	16	19	23	27	31	36	41	46	51	56	61	66	70	75	78	81

Note: This data was collected in GoldBio labs using GoldBio reagents and calculated using 100 ml volumes. All reagent volumes recorded above were adjusted accordingly to create this protocol.

Tris pKa at 25°C: 8.06  
Tris pH range: 7.0 – 9.0  
d(pKa)/dT value: -0.028

## Additional Resources

### [Biological Buffers Resource Page](#)

- Tools
- Protocols
- Articles
- Preparation Instructions
- Reference Charts



### [Stock Solution Calculator](#)

This calculator allows you estimate how much volume you'll get from a product pack size or weighed mass of powder. Simply select the product solution you're interested in, enter the mass and change units accordingly to get your volume results. This calculator also includes valuable product information including a link to the product, the stock solution recipe and basic information such as catalog number and molecular weight.



### [Helpful Articles](#)

1. [What is a biological buffer and how to choose the best buffer for your experiment](#)
2. [Why is my lysis buffer not working: How to resolve 9 lysis buffer issues that might be holding you back](#)
3. [How to prepare your most frequently used buffers](#)
4. [How to pronounce biological buffers, by GoldBio](#)



## GoldBio Buffers by Category

**Table 2. Biological Buffers**

Buffer	Catalog Number
<u>ACES</u>	A-010
<u>ADA</u>	A-780
<u>BES</u>	B-780
<u>Bicine</u>	B-785
<u>Bis-Tris</u>	B-020
<u>Boric Acid</u>	B-030
<u>CAPS</u>	C-040
<u>CHES</u>	C-870
<u>HEPES, Free Acid</u>	H-400
<u>HEPES, Sodium Salt</u>	H-401
<u>HEPPSO</u>	H-480
<u>MES, Free Acid</u>	M-095
<u>MES Free Acid Monohydrate, Ultra Pure</u>	M-090
<u>MES, Sodium Salt</u>	M-091
<u>MOPS Free Acid, Ultra Pure</u>	M-790
<u>MOPS, Sodium Salt</u>	M-791
<u>MOPSO, Free Acid</u>	M-795
<u>MOPSO, Sodium Salt</u>	M-799
<u>PBS (Phosphate Buffered Saline) Tablets</u>	P-271
<u>PIPES, Sodium Salt</u>	P-280
<u>TAPS</u>	T-780
<u>TES</u>	T-785
<u>Tricine</u>	T-870
<u>Tris Acetate</u>	T-090
<u>Tris HCl</u>	T-095
<u>Tris (Tris Base)</u>	T-400

**Table 3. Chromatography Buffers**

Buffer	Catalog Number
<u>5X Phosphate Buffer + 3M Imidazole Solution</u>	I-905
<u>5X Phosphate Buffer, pH 8</u>	P-500
<u>His-Tag Buffer Set</u>	I-906
<u>Imidazole - 3M Recrystallized Solution, pH 6.0</u>	I-901
<u>Imidazole, Recrystallized</u>	I-902

**Table 4. Cell Culture Buffers**

Buffer	Catalog Number
<u>ACES</u>	A-010
<u>BES</u>	B-780
<u>Bicine</u>	B-785
<u>HEPES, Free Acid</u>	H-400
<u>HEPES, Sodium Salt</u>	H-401
<u>MES Free Acid Monohydrate, Ultra Pure</u>	M-090
<u>MOPS Free Acid, Ultra Pure</u>	M-790
<u>MOPSO, Free Acid</u>	M-795
<u>PBS (Phosphate Buffered Saline) Tablets</u>	P-271
<u>PIPES, Sodium Salt</u>	P-280
<u>TAPS</u>	T-780
<u>Tricine</u>	T-870

**Table 5. Electrophoresis Buffers**

Buffer	Catalog Number
<u>Bis-Tris</u>	B-020
<u>Boric Acid</u>	B-030
<u>Glycine, sodium salt</u>	G-450
<u>Tricine</u>	T-870
<u>Tris Acetate</u>	T-090
<u>Tris HCl</u>	T-095
<u>Tris (Tris Base)</u>	T-400

**Table 6. Lysis Buffers**

Buffer	Catalog Number
<u>Bacterial Cell Lysis Buffer</u>	GB-176, GB-177
<u>Mammalian Cell Lysis Buffer</u>	GB-180
<u>Tissue Cell Lysis Buffer</u>	GB-181
<u>Yeast Lysis Buffer Kit plus Zymolyase</u>	GB-178
<u>Yeast Spheroplast Lysis Buffer</u>	GB-179

## Biological Buffer Selection Guide

Buffer	Catalog #	pH Range	pKa (25°C)	Applications	Metal Binding	Comments
• ACES	<a href="#">A-010</a>	6.1 – 6.8	6.8	Cell culture media, agarose and polyacrylamide electrophoresis, isoelectric focusing of proteins, x-ray crystallography, and yeast and bacterial cell studies.	Binds $\text{Cu}^{2+}$ and $\text{Mg}^{2+}$ .	Significant absorption of UV light at 230 nm. Inhibits GABA receptor binding.
• ADA	<a href="#">A-780</a>	6.0 – 7.2	6.6	Protein crystallization, electrophoresis, differential scanning calorimetry and metal decontamination in soil.	Strong binding to $\text{Mn}^{2+}$ , $\text{Co}^{2+}$ , $\text{Ni}^{2+}$ , $\text{Zn}^{2+}$ , $\text{Cd}^{2+}$ , $\text{Pb}^{2+}$ and $\text{Cu}^{2+}$ .	Absorbs UV light between 0.1 and 260 nm. Not suitable for BCA assay.
• BES	<a href="#">B-780</a>	6.4 – 7.8	7.1	Transfection, cell culture and protein quantification. Chromatography (gel filtration and cation exchange). Bacterial endotoxin studies.	Binds $\text{Cu}^{2+}$ and $\text{Co}^{2+}$ weakly.	Suitable for BCA assay. Binds DNA and interferes with restriction enzymes.
• Bicine	<a href="#">B-785</a>	7.6 – 9.0	8.3	Tissue culture, phosphorylation, protein crystallization, enzymatic reactions, a multiphasic buffer for SDS-PAGE. Cation exchange chromatography and NMR spectroscopy.	Strong binding to $\text{Mg}^{2+}$ , $\text{Ca}^{2+}$ , $\text{Co}^{2+}$ , $\text{Fe}^{3+}$ and $\text{Cu}^{2+}$ Weak binding to $\text{Mn}^{2+}$ .	Zwitterionic buffer. Not suitable for BCA assay.
• Bis-Tris	<a href="#">B-020</a>	5.8 – 7.2	6.5	Electrophoresis (sample, gel and running buffer), anion exchange chromatography, protein purification, NMR spectroscopy and x-ray crystallography.	Binds $\text{Cu}^{2+}$ and $\text{Pb}^{2+}$ . Weak binding with other metals.	Zwitterionic buffer. Substitute for cacodylate buffer. Interacts with human liver FAB and affects protein dynamics. Not suitable for BCA assay.
• Boric Acid	<a href="#">B-030</a>	8.2 – 10.2	9.2	Capillary and gel electrophoresis (protein, nucleic acid and secondary metabolite separation). Plant micronutrient.		Binds carbohydrates and pyridine nucleotides, such as $\text{NAD}^+$ .

## Biological Buffer Selection Guide (part 2)

Buffer	Catalog #	pH Range	pKa (25°C)	Applications	Metal Binding	Comments
CAPS	<a href="#">C-040</a>	9.7 - 11.1	10.4	SDS-PAGE, protein sequencing and quantification and enzymatic assays. Cation exchange chromatography and protein crystallization.	Negligible metal ion binding	Zwitterionic buffer. Suitable for BCA assay.
CHES	<a href="#">C-870</a>	8.6 - 10.0	9.5	Electrophoresis and enzymatic assays. Electron paramagnetic resonance (EPR) spectroscopy, kinetic measurements, yeast dextrose medium.	Negligible metal ion binding	Zwitterionic buffer. High affinity for iodoacetate binding site of liver alcohol dehydrogenase.
HEPES Free Acid HEPES Sodium Salt	<a href="#">H-400</a> <a href="#">H-401</a>	6.8 - 8.2	7.5	Cell culture, electroporation, qPCR, restriction enzymes, atomic force microscopy and electrophoresis. Protein quantification. Cation exchange chromatography.	No binding	Zwitterionic buffer. May form radicals and not suitable for redox reactions. Binds DNA and affects restriction enzyme function.
HEPPSO	<a href="#">H-480</a>	7.1 - 8.5	7.8	Protein quantification. Toxicology studies. Isoelectric focusing and enzymatic studies.	Weak binding with Cu <sup>2+</sup>	May form radicals and not suitable for redox reactions. Suitable for BCA assay.
MES Free Acid MES Free Acid Monohydrate, Ultra Pure MES Sodium Salt	<a href="#">M-095</a> <a href="#">M-090</a> <a href="#">M-091</a>	5.5 - 6.7	6.1	Cell and plant culture media, protein electrophoresis, chromatography, electrochromatography. Toxicology studies in yeast (metal toxicity and removal). Fluorescence microscopy.	Negligible binding	Zwitterionic buffer. Substitute for cacodylate, citrate and malate buffers. Inhibits connexin channels. Binds human liver FAB protein.
MOPS Free Acid MOPS Sodium Salt	<a href="#">M-790</a> <a href="#">M-791</a>	6.5 - 7.9	7.1	Cell culture, nucleic acid/ protein electrophoresis, and protein purification and quantification. Toxicology studies in yeast (metal toxicity).	Negligible binding	Zwitterionic buffer. Filter sterilization required. Suitable for BCA assay.

## Biological Buffer Selection Guide (part 3)

Buffer	Catalog #	pH Range	pKa (25°C)	Applications	Metal Binding	Comments
MOPSO Free Acid MOPSO Sodium Salt	<a href="#">M-795</a> <a href="#">M-799</a>	6.2 - 7.6	6.9	Cell culture media, capillary electrophoresis and quantification. Fluorescence spectroscopy, spectrophotometry and isothermal titration calorimetry. Electrochromatography.	Binds Fe <sup>3+</sup> .	Zwitterionic buffer. PTFE-coated or plastic coated utensils recommended for its use.
PIPES Free Acid PIPES Sodium Salt	<a href="#">P-281</a> <a href="#">P-280</a>	6.1 - 7.5	6.8	Cell culture, protein crystallization, electrophoresis, isoelectric focusing. Chromatography and protein quantification. Electron microscopy.	Negligible binding	Zwitterionic buffer. May form radicals and not suitable for redox reactions. Suitable for BCA assay. Substitute for cacodylate.
TES	<a href="#">T-785</a>	6.8 - 8.2	7.4	DNA precipitation and extraction. Gel filtration and chromatography. Enzyme assays. Protein quantification.	Binds strongly to Cr <sup>3+</sup> and Fe <sup>3+</sup> , and binds weakly to Co <sup>2+</sup> , Ni <sup>2+</sup> , Cu <sup>2+</sup> and Zn <sup>2+</sup>	Zwitterionic buffer. Suitable for BCA assay.
Tricine	<a href="#">T-870</a>	7.4 - 8.8	8.1	SDS-PAGE, capillary zone electrophoresis, HPLC and ion exchange chromatography, firefly luciferase assays. Lowry protein assay upon addition of	Binds Cu <sup>2+</sup> strongly, Ca <sup>2+</sup> and Mg <sup>2+</sup> moderately and weakly to Mn <sup>2+</sup>	Substitute for barbital. Photooxidized by flavins and flavoproteins.
Tris (Tris Base) Tris-HCl	<a href="#">T-400</a> <a href="#">T-095</a>	7.0 - 9.0	8.1	Electrophoresis, SDS-PAGE Laemmli buffer and other running/loading buffers with glycine, and SDS. Anion exchange chromatography.	Strong binding to Fe <sup>3+</sup> , Cr <sup>3+</sup> , Ni <sup>2+</sup> Co <sup>2+</sup> , Cu <sup>2+</sup>	Reacts with DEPC, aldehydes, common metals, various enzymes and DNA. Not suitable for BCA assay. Not suitable for most cell culture work.
TAPS	<a href="#">T-780</a>	7.7 - 9.1	8.4	DNA analysis, protein purification (ion exchange chromatography, planar chromatography), enzymatic assay (aminopeptidase activity), cell culture media (dinoflagellates).	Binds strongly to Cr <sup>3+</sup> , Fe <sup>3+</sup> and Cu <sup>2+</sup>	Inhibits connexin channels.



## Guide: Commonly used Buffers

Buffer	pH Range	Applications	Comments
PBS (Phosphate Buffered Saline)	7.3 - 7.5	Cell culture, immunoassays, sample dilution and protein purification.	Isotonic and nontoxic buffer. Mimics pH, osmolarity and ion concentration of the human body. pKa 7.4. <a href="#">GoldBio Catalog # P-271</a> .
Cacodylate	5.0 - 7.4	Electron microscopy and histology.	Substitute for Sorensen's buffer. Toxic. At low pH, oxidizes thiol groups and can inactivate enzymes that require reduced thiols. pKa 6.27.
Stripping buffer		Western blotting.	Components: Tris-HCl 1M, 20% SDS and $\beta$ -mercaptoethanol. Prepare fresh prior to use. If antibody recovery from the membrane is desired, avoid SDS.
Transfer buffer		Western blotting.	Components: Tris-base 25mM, glycine 192mM, and 20% methanol.
Running buffer	8.3	Western blotting.	Components: Tris-base 25mM, glycine 192mM, and 0.1% SDS. Can be stored at room temperature. No pH adjustment is required.
Krebs-Henseleit bicarbonate buffer	7.4	Perfusion and isolation of muscle tissue.	Components: NaCl 118mM, KCl 4.7mM, MgSO <sub>4</sub> 1.2mM, CaCl <sub>2</sub> 2.5mM and NaHCO <sub>3</sub> 25mM. Equilibrate with 95% O <sub>2</sub> and 5% CO <sub>2</sub> . Add CaCl <sub>2</sub> after all other compounds are dissolved. Add glucose for cell maintenance. Use immediately after preparation.
TE (Tris-EDTA buffer)	7.9 - 8.1	Storage of nucleic acids.	Components: Tris-Cl 10mM pH 7.4-8.0, EDTA 1mM pH 8.0. EDTA might affect downstream applications. Can be stored at room temperature for up to 6 months.

## Guide: Commonly used Buffers (part 2)

Buffer	pH Range	Applications	Comments
TAE (Tris-acetate-EDTA buffer)	8.0 - 8.3	Agarose electrophoresis for the separation of nucleic acids (DNA and RNA).	Components: Tris base 40mM, acetic acid 20mM and 1mM EDTA. TAE has a lower buffering capacity than TBE. Can be stored at room temperature. <a href="#">Stock solution protocol.</a>
TTE (Tris-TAPS-EDTA buffer)	8.3	Electrophoresis.	Components: 500mM Tris base, 500mM TAPS and 10mM EDTA. Can be stored at room temperature for up to 4 weeks or at 4°C for long term storage. <a href="#">Stock solution protocol.</a>
TBE (Tris-borate-EDTA buffer)	8.3	Polyacrylamide gel electrophoresis for the separation of nucleic acids (DNA and RNA). blotting.	Components: Tris base 89mM, boric acid 89mM and EDTA 2mM. Used with denaturing and non-denaturing gels. Borate strongly inhibits enzymatic activity. <a href="#">Stock solution protocol.</a>
TBST (Tris buffered saline with Tween® 20)	7.4 - 7.6	Western blotting and ELISA.	Components: Tris HCl 20mM pH 7.5, NaCl 150mM and 0.05-0.1% Tween® 20.
Tris-Glycine	8.3	Western blotting, SDS-PAGE, ion-exchange chromatography.	Components: Tris-base 25mM, glycine 250mM and 0.1% SDS.
Imidazole-HCl	6.2 - 7.8	Affinity chromatography and SDS-PAGE.	Binds various metals. Unstable.

# The Ultimate Buffer Selection Guide

-arranged by pH

pH Range: 5 | 6 | 7 | 8 | 9 | 10 | 11

Buffer:

**MES (M-090)**

**Bis-Tris (B-020)**

**ADA (A-780)**

**ACES (A-010)**

**PIPES (P-281)**

**MOPSO (M-795)**

**BES (B-780)**

**MOPS (M-790)**

**HEPES (H-400)**

**TES (T-785)**

**Tris (T-400)**

**HEPPSO (H-480)**

**PBS (P-271)**

**Tricine (T-870)**


**Bicine (B-785)**

**TAPS (T-780)**

**Boric Acid (B-031)**

**CHES (C-870)**

**CAPS (C-040)**

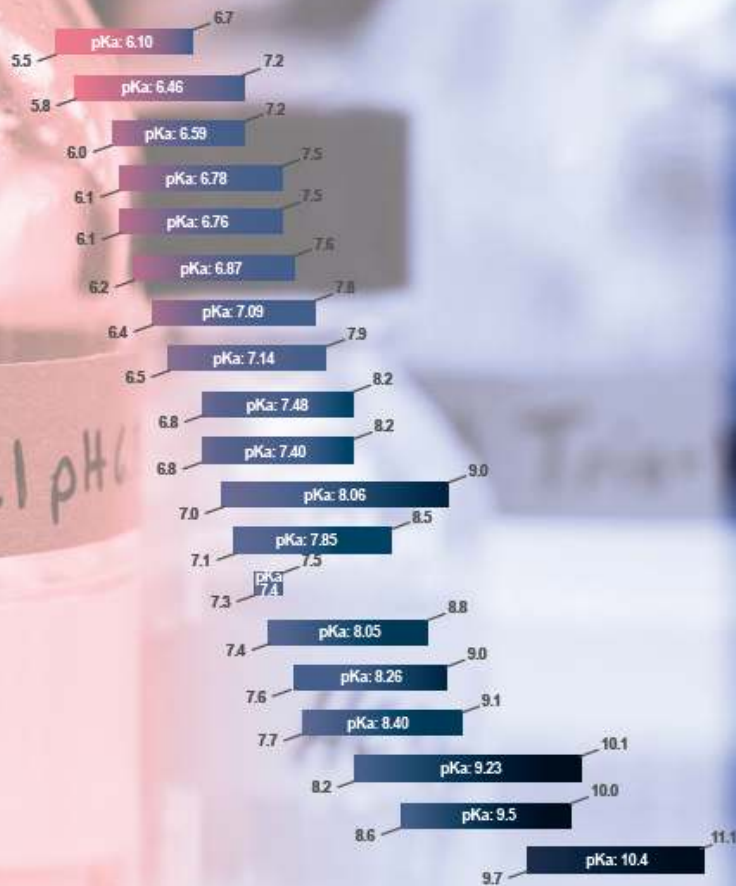

## References

- Ahmad, A., & Ghasemi, J. (2007). New buffers to improve the quantitative real-time polymerase chain reaction. *Bioscience, Biotechnology, and Biochemistry*, 71(8), 1970-1978.
- Ahmedna, M., Clarke, S. J., Rao, R. M., Marshall, W. E., & Johns, M. M. (1997). Use of filtration and buffers in raw sugar colour measurements. *Journal of the Science of Food and Agriculture*, 75(1), 109-116.
- Alonso, A. D. C., Zaidi, T., Novak, M., Grundke-Iqbal, I., & Iqbal, K. (2001). Hyperphosphorylation induces self-assembly of  $\tau$  into tangles of paired helical filaments/straight filaments. *Proceedings of the National Academy of Sciences*, 98(12), 6923-6928.
- Anderson, M. P., & Gronwald, J. W. (1991). Atrazine resistance in a velvetleaf (*Abutilon theophrasti*) biotype due to enhanced glutathione S-transferase activity. *Plant Physiology*, 96(1), 104-109.
- Antonetti, D. A., Barber, A. J., Hollinger, L. A., Wolpert, E. B., & Gardner, T. W. (1999). Vascular endothelial growth factor induces rapid phosphorylation of tight junction proteins occludin and zonula occluden 1 A potential mechanism for vascular permeability in diabetic retinopathy and tumors. *Journal of Biological Chemistry*, 274(33), 23463-23467.
- Aronoff, S., Chen, T. C., & Cheveldayoff, M. (1975). Complexation of D-glucose with borate. *Carbohydrate Research*, 40(2), 299-309.
- Baker, C. J., Mock, N. M., Roberts, D. P., Deahl, K. L., Hapeman, C. J., Schmidt, W. F., & Kochansky, J. (2007). Interference by Mes [2-(4-morpholino) ethanesulfonic acid] and related buffers with phenolic oxidation by peroxidase. *Free Radical Biology and Medicine*, 43(9), 1322-1327.
- Baker, C. J., Mock, N. M., Roberts, D. P., Deahl, K. L., Hapeman, C. J., Schmidt, W. F., & Kochansky, J. (2007). Interference by Mes [2-(4-morpholino) ethanesulfonic acid] and related buffers with phenolic oxidation by peroxidase. *Free Radical Biology and Medicine*, 43(9), 1322-1327.
- Basto, S., Dorca-Fornell, C., Thompson, K., & Rees, M. (2013). Effect of pH buffer solutions on seed germination of *Hypericum pulchrum*, *Campanula rotundifolia* and *Scabiosa columbaria*. *Seed Science and Technology*, 41(2), 298-302.
- Baur, P. S., & Stacey, T. R. (1977). The use of PIPES buffer in the fixation of mammalian and marine tissues for electron microscopy. *Journal of Microscopy*, 109(3), 315-327.
- Behr, B. R., Stratton, C. J., Foote, W. D., Knutzen, V., & Sher, G. (1990). In vitro fertilization (IVF) of mouse ova in HEPES-buffered culture media. *Journal of in Vitro Fertilization and Embryo Transfer*, 7(1), 9-15.
- Bevans, C. G., & Harris, A. L. (1999). Regulation of Connexin Channels by pH Direct Action of the Protonated Form of Taurine and other Aminosulfonates. *Journal of Biological Chemistry*, 274(6), 3711-3719.
- Blaber, S. I., Culajay, J. F., Khurana, A., & Blaber, M. (1999). Reversible thermal denaturation of human FGF-1 induced by low concentrations of guanidine hydrochloride. *Biophysical Journal*, 77(1), 470-477.

- Boughtflower, R. J., Underwood, T., & Paterson, C. J. (1995). Capillary electrochromatography—Some important considerations in the preparation of packed capillaries and the choice of mobile phase buffers. *Chromatographia*, *40*(5-6), 329-335.
- Bouwmeester, H. J., Wallaart, T. E., Janssen, M. H., van Loo, B., Jansen, B. J., Posthumus, M. A., ... & Franssen, M. C. (1999). Amorpha-4, 11-diene synthase catalyses the first probable step in artemisinin biosynthesis. *Phytochemistry*, *52*(5), 843-854.
- Brignac Jr, P. J., & Mo, C. (1975). Formation constants and metal-to-ligand ratios for tris (hydroxymethyl) aminomethane-metal complexes. *Analytical Chemistry*, *47*(8), 1465-1466.
- Brody, J. R., & Kern, S. E. (2004). History and principles of conductive media for standard DNA electrophoresis. *Analytical Biochemistry*, *333*(1), 1-13.
- Brown, R. E., Jarvis, K. L., & Hyland, K. J. (1989). Protein measurement using bicinchoninic acid: elimination of interfering substances. *Analytical Biochemistry*, *180*(1), 136-139.
- Budowle, B. (1987). Improved separation of the common transferrin variants in gels containing pH 5–7 Ampholines and HEPES. *Electrophoresis*, *8*(4), 210-212.
- Celie, P. H., Klaassen, R. V., van Rossum-Fikkert, S. E., van Elk, R., van Nierop, P., Smit, A. B., & Sixma, T. K. (2005). Crystal structure of acetylcholine-binding protein from *Bulinus truncatus* reveals the conserved structural scaffold and sites of variation in nicotinic acetylcholine receptors. *Journal of Biological Chemistry*, *280*(28), 26457-26466.
- Chen, C., & Okayama, H. (1987). High-efficiency transformation of mammalian cells by plasmid DNA. *Molecular and Cellular Biology*, *7*(8), 2745-2752.
- Chen, T. C., & Hong, A. (1995). Chelating extraction of lead and copper from an authentic contaminated soil using N-(2-acetamido) iminodiacetic acid and S-carboxymethyl-L-cysteine. *Journal of Hazardous Materials*, *41*(2), 147-160.
- Clayton, L., Quinlan, R. A., Roobol, A., Pogson, C. I., & Gull, K. (1980). A comparison of tubulins from mammalian brain and *Physarum polycephalum* using SDS—polyacrylamide gel electrophoresis and peptide mapping. *FEBS Letters*, *115*(2), 301-305.
- Cockcroft, S., Thomas, G. L., Fensome, A., Geny, B., Cunningham, E., Gout, I., ... & Hsuan, J. J. (1994). Phospholipase D: a downstream effector of ARF in granulocytes. *Science*, *263*(5146), 523-526.
- Colby, S. M., Alonso, W. R., Katahira, E. J., McGarvey, D. J., & Croteau, R. (1993). 4S-limonene synthase from the oil glands of spearmint (*Mentha spicata*). cDNA isolation, characterization, and bacterial expression of the catalytically active monoterpene cyclase. *Journal of Biological Chemistry*, *268*(31), 23016-23024.
- Collins, C. A., Fry, F. H., Holme, A. L., Yiakouvaki, A., Al-Qenaei, A., Pourzand, C., & Jacob, C. (2005). Towards multifunctional antioxidants: synthesis, electrochemistry, in vitro and cell culture evaluation of compounds with ligand/catalytic properties. *Organic & Biomolecular Chemistry*, *3*(8), 1541-1546.
- Comai, L. U. C. A., & Kosuge, T. S. U. N. E. (1980). Involvement of plasmid deoxyribonucleic acid in indoleacetic acid synthesis in *Pseudomonas savastanoi*. *Journal of Bacteriology*, *143*(2), 950-957.

- Cook, T. A., Urrutia, R., & McNiven, M. A. (1994). Identification of dynamin 2, an isoform ubiquitously expressed in rat tissues. *Proceedings of the National Academy of Sciences*, 91(2), 644-648.
- Cunliffe, J. M., Baryla, N. E., & Lucy, C. A. (2002). Phospholipid bilayer coatings for the separation of proteins in capillary electrophoresis. *Analytical Chemistry*, 74(4), 776-783.
- Dahl, C., Schulte, A., & Shin, D. H. (2007). Cloning, expression, purification, crystallization and preliminary X-ray diffraction analysis of DsrEFH from *Allochrocatium vinosum*. *Acta Crystallographica Section F: Structural Biology and Crystallization Communications*, 63(10), 890-892.
- de Souza, M. L., Sadowsky, M. J., & Wackett, L. P. (1996). Atrazine chlorohydrolase from *Pseudomonas* sp. strain ADP: gene sequence, enzyme purification, and protein characterization. *Journal of Bacteriology*, 178(16), 4894-4900.
- Deutscher, M. P. (Ed.). (1990). *Guide to Protein Purification* (Vol. 182). Gulf Professional Publishing.
- Donarski, W. J., Dumas, D. P., Heitmeyer, D. P., Lewis, V. E., & Raushel, F. M. (1989). Structure-activity relationships in the hydrolysis of substrates by the phosphotriesterase from *Pseudomonas diminuta*. *Biochemistry*, 28(11), 4650-4655.
- Donoghue, M., Hsieh, F., Baronas, E., Godbout, K., Gosselin, M., Stagliano, N., ... & Acton, S. (2000). A novel angiotensin-converting enzyme-related carboxypeptidase (ACE2) converts angiotensin I to angiotensin 1-9. *Circulation Research*, 87(5), e1-e9.
- Downs, S. M., & Mastropolo, A. M. (1997). Culture conditions affect meiotic regulation in cumulus cell-enclosed mouse oocytes. *Molecular Reproduction and Development*, 46(4), 551-566.
- Edelstein, P. H., & Edelstein, M. A. (1993). Comparison of three buffers used in the formulation of buffered charcoal yeast extract medium. *Journal of Clinical Microbiology*, 31(12), 3329-3330.
- El-Harakany, A. A., Halim, F. A., & Barakat, A. O. (1984). Dissociation constants and related thermodynamic quantities of the protonated acid form of tris-(hydroxymethyl)-aminomethane in mixtures of 2-methoxyethanol and water at different temperatures. *Journal of Electroanalytical Chemistry and Interfacial Electrochemistry*, 162(1), 285-305.
- Ferguson, W. J., Braunschweiger, K. I., Braunschweiger, W. R., Smith, J. R., McCormick, J. J., Wasmann, C. C., ... & Good, N. E. (1980). Hydrogen ion buffers for biological research. *Analytical Biochemistry*, 104(2), 300-310.
- Ferreira, C. M., Pinto, I. S., Soares, E. V., & Soares, H. M. (2015). (Un) suitability of the use of pH buffers in biological, biochemical and environmental studies and their interaction with metal ions—a review. *RSC Advances*, 5(39), 30989-31003.
- Fromm, M., Taylor, L. P., & Walbot, V. (1985). Expression of genes transferred into monocot and dicot plant cells by electroporation. *Proceedings of the National Academy of Sciences*, 82(17), 5824-5828.
- Funtenberger, S., Dumay, E., & Cheftel, J. C. (1995). Pressure-induced aggregation of  $\beta$ -lactoglobulin in pH 7.0 buffers. *LWT-Food Science and Technology*, 28(4), 410-418.



- Galant, A., Arkus, K. A., Zubieta, C., Cahoon, R. E., & Jez, J. M. (2009). Structural basis for evolution of product diversity in soybean glutathione biosynthesis. *The Plant Cell*, 21(11), 3450-3458.
- Gardner, R. S. (1969). The use of tricine buffer in animal tissue cultures. *The Journal of Cell Biology*, 42(1), 320-321.
- Good, N. E., & Izawa, S. (1972). Hydrogen ion buffers. *Methods in Enzymology*, 24, 53.
- Good, N. E., Winget, G. D., Winter, W., Connolly, T. N., Izawa, S., & Singh, R. M. (1966). Hydrogen ion buffers for biological research\*. *Biochemistry*, 5(2), 467-477.
- Grady, J. K., Chasteen, N. D., & Harris, D. C. (1988). Radicals from "Good's" buffers. *Analytical Biochemistry*, 173(1), 111-115.
- Grande, H. J., & Van Der Ploeg, K. R. (1978). Tricine radicals as formed in the presence of peroxide producing enzymes. *FEBS Letters*, 95(2), 352-356.
- Greenway, H., & Osmond, C. B. (1972). Salt responses of enzymes from species differing in salt tolerance. *Plant Physiology*, 49(2), 256-259.
- Habib, A., & Tabata, M. (2004). Oxidative DNA damage induced by HEPES (2-[4-(2-hydroxyethyl)-1-piperazinyl] ethanesulfonic acid) buffer in the presence of Au (III). *Journal of Inorganic Biochemistry*, 98(11), 1696-1702.
- Hammadeh, M. E., Greiner, S., Rosenbaum, P., & Schmidt, W. (2001). Comparison between human sperm preservation medium and TEST-Yolk buffer on protecting chromatin and morphology integrity of human spermatozoa in fertile and subfertile men after freeze-thawing procedure. *Journal of Andrology*, 22(6), 1012-1018.
- Hamoudová, R., & Pospíšilová, M. (2006). Determination of ibuprofen and flurbiprofen in pharmaceuticals by capillary zone electrophoresis. *Journal of Pharmaceutical and Biomedical Analysis*, 41(4), 1463-1467.
- Hicks, M., & Gebicki, J. M. (1986). Rate constants for reaction of hydroxyl radicals with Tris, Tricine and HEPES buffers. *FEBS Letters*, 199(1), 92-94.
- Hoskins, L. C., Agustines, M., McKee, W. B., Boulding, E. T., Kriaris, M., & Niedermeyer, G. (1985). Mucin degradation in human colon ecosystems. Isolation and properties of fecal strains that degrade ABH blood group antigens and oligosaccharides from mucin glycoproteins. *Journal of Clinical Investigation*, 75(3), 944.
- Huisman, T. H., & Dozy, A. M. (1965). Studies on the heterogeneity of hemoglobin: IX. The use of tris (hydroxymethyl) aminomethane—HCl buffers in the anion-exchange chromatography of hemoglobins. *Journal of Chromatography A*, 19, 160-169.
- Hulett, F. M., Bookstein, C., & Jensen, K. (1990). Evidence for two structural genes for alkaline phosphatase in *Bacillus subtilis*. *Journal of Bacteriology*, 172(2), 735-740.
- Innis, M. A., Gelfand, D. H., & Sninsky, J. J. (Eds.). (1999). *PCR Applications: Protocols for Functional Genomics*. Academic press.

- Itagaki, A., & Kimura, G. (1974). TES and HEPES buffers in mammalian cell cultures and viral studies: problem of carbon dioxide requirement. *Experimental Cell Research*, 83(2), 351-361.
- Jennings, C., West, J., Waine, C., Craik, D., & Anderson, M. (2001). Biosynthesis and insecticidal properties of plant cyclotides: the cyclic knotted proteins from *Oldenlandia affinis*. *Proceedings of the National Academy of Sciences*, 98(19), 10614-10619.
- Jensen, P. R., & Hammer, K. (1993). Minimal requirements for exponential growth of *Lactococcus lactis*. *Applied and Environmental Microbiology*, 59(12), 4363-4366.
- Johnson, S. L., & Smith, K. W. (1976). The interaction of borate and sulfite with pyridine nucleotides. *Biochemistry*, 15(3), 553-559.
- Kallweit, U., Börnsen, K. O., Kresbach, G. M., & Widmer, H. M. (1996). Matrix compatible buffers for analysis of proteins with matrix-assisted laser desorption/ionization mass spectrometry. *Rapid Communications in Mass Spectrometry*, 10(7), 845-849.
- Kaneta, T., Tanaka, S., & Yoshida, H. (1991). Improvement of resolution in the capillary electrophoretic separation of catecholamines by complex formation with boric acid and control of electroosmosis with a cationic surfactant. *Journal of Chromatography A*, 538(2), 385-391.
- Kanyo, Z. F., Chen, C. Y., Daghigh, F., Ash, D. E., & Christianson, D. W. (1992). Crystallization and oligomeric structure of rat liver arginase. *Journal of Molecular Biology*, 224(4), 1175-1177.
- Kashino, Y., Koike, H., & Satoh, K. (2001). An improved sodium dodecyl sulfate-polyacrylamide gel electrophoresis system for the analysis of membrane protein complexes. *Electrophoresis*, 22(6), 1004-1007.
- Kaushal, V., & Barnes, L. D. (1986). Effect of zwitterionic buffers on measurement of small masses of protein with bicinchoninic acid. *Analytical Biochemistry*, 157(2), 291-294.
- Keating, K. I., Caffrey, P. B., & Dagbusan, B. C. (1996). Buffers in daphnid culture and bioassay. *Environmental Toxicology and Chemistry*, 15(3), 348-352.
- Kelly, A. E., Ou, H. D., Withers, R., & Dötsch, V. (2002). Low-conductivity buffers for high-sensitivity NMR measurements. *Journal of the American Chemical Society*, 124(40), 12013-12019.
- Kingston, R. E. (1991). Introduction of DNA into mammalian cells. *Current Protocols in Molecular Biology*, 9-0.
- Kirsch, M., Lomonosova, E. E., Korth, H. G., Sustmann, R., & de Groot, H. (1998). Hydrogen Peroxide Formation by Reaction of Peroxynitrite with HEPES and Related Tertiary Amines Implications for a General Mechanism. *Journal of Biological Chemistry*, 273(21), 12716-12724.
- Klampfl, C. W., Katzmayer, M. U., & Buchberger, W. (1998). Separation of inorganic and organic anions by capillary zone electrophoresis with simultaneous indirect UV and conductivity detection. *Electrophoresis*, 19(14), 2459-2464.
- Kohoutova, J., Kutá Smatanová, I., Brynda, J., Lapkouski, M., Revuelta, J. L., Arellano, J. B., & Ettrich, R. (2009). Crystallization and preliminary crystallographic characterization of the extrinsic PsbP protein of photosystem II from *Spinacia oleracea*. *Acta Crystallographica Section F: Structural Biology and Crystallization Communications*, 65(2), 111-115.



- Koort, J. M., Lukinmaa, S., Rantala, M., Unkila, E., & Siitonen, A. (2002). Technical improvement to prevent DNA degradation of enteric pathogens in pulsed-field gel electrophoresis. *Journal of Clinical Microbiology*, 40(9), 3497-3498.
- Lavigne, V., Pons, A., & Dubourdieu, D. (2007). Assay of glutathione in must and wines using capillary electrophoresis and laser-induced fluorescence detection: changes in concentration in dry white wines during alcoholic fermentation and aging. *Journal of Chromatography A*, 1139(1), 130-135.
- Littlechild, J., Turner, N., Hobbs, G., Lilly, M., Rawas, A., & Watson, H. (1995). Crystallization and preliminary X-ray crystallographic data with *Escherichia coli* transketolase. *Acta Crystallographica Section D: Biological Crystallography*, 51(6), 1074-1076.
- Liu, Q., Li, X., & Sommer, S. S. (1999). pK-matched running buffers for gel electrophoresis. *Analytical Biochemistry*, 270(1), 112-122.
- Long, D., & Yang, D. (2009). Buffer interference with protein dynamics: a case study on human liver fatty acid binding protein. *Biophysical Journal*, 96(4), 1482-1488.
- Lu, Z., Breidt, F., Fleming, H. P., Altermann, E., & Klaenhammer, T. R. (2003). Isolation and characterization of a *Lactobacillus plantarum* bacteriophage, ΦJL-1, from a cucumber fermentation. *International Journal of Food Microbiology*, 84(2), 225-235.
- Lyubchenko, Y. L., Shlyakhtenko, L. S., & Ando, T. (2011). Imaging of nucleic acids with atomic force microscopy. *Methods*, 54(2), 274-283.
- Mahadevan, M. M., Fleetham, J., Church, R. B., & Taylor, P. J. (1986). Growth of mouse embryos in bicarbonate media buffered by carbon dioxide, HEPES, or phosphate. *Journal of In Vitro Fertilization and Embryo Transfer*, 3(5), 304-308.
- Marina, M. L., Benito, I., Diez-Masa, J. C., & González, M. J. (1996). Chiral separation of polychlorinated biphenyls by micellar electrokinetic chromatography with  $\gamma$ -cyclodextrin as modifier in the separation buffer. *Chromatographia*, 42(5-6), 269-272.
- Michels, D. A., Hu, S., Schoenherr, R. M., Eggertson, M. J., & Dovichi, N. J. (2002). Fully automated two-dimensional capillary electrophoresis for high sensitivity protein analysis. *Molecular & Cellular Proteomics*, 1(1), 69-74.
- Mohamed, M. M., Abd-Alla, E. M., & El-Badawy, A. E. S. (2007). Dimethyltin (IV) complexes with zwitterionic buffers (MES and MOPS). *Journal of Organometallic Chemistry*, 692(8), 1735-1747.
- Monthony, J. F., Wallace, E. G., & Allen, D. M. (1978). A non-barbital buffer for immunoelectrophoresis and zone electrophoresis in agarose gels. *Clinical Chemistry*, 24(10), 1825-1827.
- Moore, S. A., Kingston, R. L., Loomes, K. M., Hernell, O., Bläckberg, L., Baker, H. M., & Baker, E. N. (2001). The structure of truncated recombinant human bile salt-stimulated lipase reveals bile salt-independent conformational flexibility at the active-site loop and provides insights into heparin binding. *Journal of Molecular Biology*, 312(3), 511-523.

- Morgan, J. A., Barney, C. S., Penn, A. H., & Shanks, J. V. (2000). Effects of buffered media upon growth and alkaloid production of *Catharanthus roseus* hairy roots. *Applied Microbiology and Biotechnology*, 53(3), 262-265.
- Mumford, R. A., Pierzchala, P. A., Strauss, A. W., & Zimmerman, M. (1981). Purification of a membrane-bound metalloendopeptidase from porcine kidney that degrades peptide hormones. *Proceedings of the National Academy of Sciences*, 78(11), 6623-6627.
- Neidhardt, F. C., Bloch, P. L., & Smith, D. F. (1974). Culture medium for enterobacteria. *Journal of Bacteriology*, 119(3), 736-747.
- Neill, S. J., Desikan, R., Clarke, A., & Hancock, J. T. (2002). Nitric oxide is a novel component of abscisic acid signaling in stomatal guard cells. *Plant Physiology*, 128(1), 13-16.
- Nijtmans, L. G., Henderson, N. S., & Holt, I. J. (2002). Blue Native electrophoresis to study mitochondrial and other protein complexes. *Methods*, 26(4), 327-334.
- Nishiwaki, T., & Hayashi, K. (2001). Purification and characterization of an aminopeptidase from the edible basidiomycete *Grifola frondosa*. *Bioscience, Biotechnology, and Biochemistry*, 65(2), 424-427.
- Nurok, D., Frost, M. C., & Chenoweth, D. M. (2000). Separation using planar chromatography with electroosmotic flow. *Journal of Chromatography A*, 903(1), 211-217.
- Omburo, G. A., Kuo, J. M., Mullins, L. S., & Raushel, F. (1992). Characterization of the zinc binding site of bacterial phosphotriesterase. *Journal of Biological Chemistry*, 267(19), 13278-13283.
- Palasz, A. T., Breña, P. B., De la Fuente, J., & Gutierrez-Adan, A. (2008). The effect of different zwitterionic buffers and PBS used for out-of-incubator procedures during standard in vitro embryo production on development, morphology and gene expression of bovine embryos. *Theriogenology*, 70(9), 1461-1470.
- Parfitt, D. E., Almehdi, A. A., & Bloksberg, L. N. (1988). Use of organic buffers in plant tissue-culture systems. *Scientia Horticulturae*, 36(3), 157-163.
- Perez-Miller, S. J., & Hurley, T. D. (2003). Coenzyme isomerization is integral to catalysis in aldehyde dehydrogenase. *Biochemistry*, 42(23), 7100-7109.
- Pérez-Ruiz, T., Martínez-Lozano, C., Sanz, A., & Bravo, E. (2002). Development and validation of a capillary electrophoretic method for the determination of amiodarone and desethylamiodarone. *Chromatographia*, 56(1-2), 63-67.
- Peterson, G. L. (1979). Review of the Folin phenol protein quantitation method of Lowry, Rosebrough, Farr and Randall. *Analytical Biochemistry*, 100(2), 201-220.
- Poole, C. A., Reilly, H. C., & Flint, M. H. (1982). The adverse effects of HEPES, TES, and BES zwitterion buffers on the ultrastructure of cultured chick embryo epiphyseal chondrocytes. *In vitro*, 18(9), 755-765.
- Prange, A., & Schaumlöffel, D. (1999). Determination of element species at trace levels using capillary electrophoresis-inductively coupled plasma sector field mass spectrometry. *Journal of Analytical Atomic Spectrometry*, 14(9), 1329-1332.

- Praveen, V. (2009, November 18). *Tris-HEPES SDS-PAGE*. Retrieved August 5, 2015, from <http://biomedicalprotocols.blogspot.com/2009/11/tris-hepes-sds-page.html>
- R Loeblich, A. (1975). A Seawater Medium for Dinoflagellates and the Nutrition of *Cachonina NIE11*. *Journal of Phycology*, 11(1), 80-86.
- Ream, P. J., Suljak, S. W., Ewing, A. G., & Han, K. A. (2003). Micellar Electrokinetic Capillary Chromatography-Electrochemical Detection for Analysis of Biogenic Amines in *Drosophila melanogaster*. *Analytical Chemistry*, 75(16), 3972-3978.
- Reichel, C., Kulovics, R., Jordan, V., Watzinger, M., & Geisendorfer, T. (2009). SDS-PAGE of recombinant and endogenous erythropoietins: benefits and limitations of the method for application in doping control. *Drug Testing and Analysis*, 1(1), 43-50.
- Richter, A. (1967). Plaque formation with N-tris (hydroxymethyl) methyl-2-aminoethanesulfonic acid (TES buffer). *Applied Microbiology*, 15(6), 1507.
- Robinson, J. D., & Davis, R. L. (1987). Buffer, pH, and ionic strength effects on the (Na<sup>++</sup> K<sup>+</sup>)-ATPase. *Biochimica et Biophysica Acta (BBA)-Protein Structure and Molecular Enzymology*, 912(3), 343-347.
- Rodriguez-Valera, F., Ruiz-Berraquero, F., & Ramos-Cormenzana, A. (1980). Short Communication Isolation of Extremely Halophilic Bacteria Able to Grow in Defined Inorganic Media with Single Carbon Sources. *Journal of General Microbiology*, 119(2), 535-538.
- Salentine, C. G. (1983). High-field boron-11 NMR of alkali borates. Aqueous polyborate equilibria. *Inorganic Chemistry*, 22(26), 3920-3924.
- Schägger, H. (2006). Tricine-SDS-PAGE. *Nature Protocols*, 1, 16-22.
- Schägger, H., & Von Jagow, G. (1987). Tricine-sodium dodecyl sulfate-polyacrylamide gel electrophoresis for the separation of proteins in the range from 1 to 100 kDa. *Analytical Biochemistry*, 166(2), 368-379.
- Schenborn, E. T., & Goiffon, V. (2000). Calcium phosphate transfection of mammalian cultured cells. *Transcription Factor Protocols* (pp. 135-145). Humana Press.
- Schirch, V., Hopkins, S., Villar, E., & Angelaccio, S. (1985). Serine hydroxymethyltransferase from *Escherichia coli*: purification and properties. *Journal of Bacteriology*, 163(1), 1-7.
- Schliwa, M., & Van Blerkom, J. (1981). Structural interaction of cytoskeletal components. *The Journal of Cell Biology*, 90(1), 222-235.
- Scopes, R. K. (2013). *Protein Purification: Principles and Practice*. Springer Science & Business Media.
- Sheng, F., Jia, X., Yep, A., Preiss, J., & Geiger, J. H. (2009). The crystal structures of the open and catalytically competent closed conformation of *Escherichia coli* glycogen synthase. *Journal of Biological Chemistry*, 284(26), 17796-17807.
- Shiraishi, H., Kataoka, M., Morita, Y., & Umemoto, J. (1993). Interactions of hydroxyl radicals with tris (hydroxymethyl) aminomethane and Good's buffers containing hydroxymethyl or hydroxyethyl residues produce formaldehyde. *Free Radical Research*, 19(5), 315-321.

- Smelt, J. P., Otten, G. D., & Bos, A. P. (2002). Modelling the effect of sublethal injury on the distribution of the lag times of individual cells of *Lactobacillus plantarum*. *International Journal of Food Microbiology*, 73(2), 207-212.
- Soltis, D. E., Haufler, C. H., Darrow, D. C., & Gastony, G. J. (1983). Starch gel electrophoresis of ferns: a compilation of grinding buffers, gel and electrode buffers, and staining schedules. *American Fern Journal*, 9-27.
- Spendlove, R. S., Crosbie, R. B., Hayes, S. F., & Keeler, R. Y. (1971). Tricine-buffered tissue culture media for control of mycoplasma contaminants. *Experimental Biology and Medicine*, 137(1), 258-263.
- Stellwagen, N. C., Bossi, A., Gelfi, C., & Righetti, P. G. (2000). DNA and buffers: are there any noninteracting, neutral pH buffers? *Analytical biochemistry*, 287(1), 167-175.
- Stoien, J. D., & Wang, R. J. (1974). Effect of near-ultraviolet and visible light on mammalian cells in culture II. Formation of toxic photoproducts in tissue culture medium by blacklight. *Proceedings of the National Academy of Sciences*, 71(10), 3961-3965.
- Susin, S. A., Lorenzo, H. K., Zamzami, N., Marzo, I., Brenner, C., Larochette, N., ... & Kroemer, G. (1999). Mitochondrial release of caspase-2 and-9 during the apoptotic process. *The Journal of Experimental Medicine*, 189(2), 381-394.
- Syvetsen, C., & McKinley-McKee, J. S. (1981). Affinity Labeling of Liver Alcohol Dehydrogenase. *European Journal of Biochemistry*, 117(1), 165-170.
- Taha, M., Gupta, B. S., Khoiroh, I., & Lee, M. J. (2011). Interactions of biological buffers with macromolecules: The ubiquitous “smart” polymer PNIPAM and the biological buffers MES, MOPS, and MOPSO. *Macromolecules*, 44(21), 8575-8589.
- Tang, P. (2005). Boric Acid Catalyzed Amide Formation from Carboxylic Acids and Amines: N-Benzyl-4-phenylbutyramide. *Organic Syntheses*, 262-272.
- Thiel, T., Liczkowski, L., & Bissen, S. T. (1998). New zwitterionic butanesulfonic acids that extend the alkaline range of four families of good buffers: Evaluation for use in biological systems. *Journal of Biochemical and Biophysical Methods*, 37(3), 117-129.
- Thomas, B. R., Vekilov, P. G., & Rosenberger, F. (1996). Heterogeneity determination and purification of commercial hen egg-white lysozyme. *Acta Crystallographica Section D: Biological Crystallography*, 52(4), 776-784.
- Trejo, G., Frausto-Reyes, C., Gama, S. C., Meas, Y., & Orozco, G. (2012). Raman Study of Benzylideneacetone on Silver. *Int. J. Electrochem. Sci*, 7, 8436-8443.
- Tunncliffe, G. and Smith, J. A. (1981), Competitive Inhibition of  $\gamma$ -Aminobutyric Acid Receptor Binding by N-2-Hydroxyethylpiperazine-N'-2-Ethanesulfonic Acid and Related Buffers. *Journal of Neurochemistry*, 36: 1122–1126.
- Ukaegbu, U. E., Kantz, A., Beaton, M., Gassner, G. T., & Rosenzweig, A. C. (2010). Structure and ligand binding properties of the epoxidase component of styrene monooxygenase. *Biochemistry*, 49(8), 1678-1688.

- Vega, C. A., Butler, R. A., Perez, B., & Torres, C. (1985). Thermodynamics of the dissociation of protonated tris (hydroxymethyl) aminomethane in 25 and 50 wt % 2-propanol from 5 to 45 degree C. *Journal of Chemical and Engineering Data*, 30(4), 376-379.
- Warrington, K. (1923). The effect of boric acid and borax on the broad bean and certain other plants. *Ann. Bot.*, 37(4), 629-672.
- Wasserman, S. A., Daub, E., Grisafi, P., Botstein, D., & Walsh, C. T. (1984). Catabolic alanine racemase from *Salmonella typhimurium*: DNA sequence, enzyme purification, and characterization. *Biochemistry*, 23(22), 5182-5187.
- Webster, J. J., Chang, J. C., Manley, E. R., Spivey, H. O., & Leach, F. R. (1980). Buffer effects on ATP analysis by firefly luciferase. *Analytical Biochemistry*, 106(1), 7-11.
- WEIDEL, L., & PRINS, G. S. (1987). Cryosurvival of human spermatozoa frozen in eight different buffer systems. *Journal of Andrology*, 8(1), 41-47.
- Weisenberg, R. C. (1972). Microtubule formation in vitro in solutions containing low calcium concentrations. *Science*, 177(4054), 1104-1105.
- Wenner, J. R., & Bloomfield, V. A. (1999). Buffer effects on EcoRV kinetics as measured by fluorescent staining and digital imaging of plasmid cleavage. *Analytical Biochemistry*, 268(2), 201-212.
- Wigler, M., Pellicer, A., Silverstein, S., Axel, R., Urlaub, G., & Chasin, L. (1979). DNA-mediated transfer of the adenine phosphoribosyltransferase locus into mammalian cells. *Proceedings of the National Academy of Sciences*, 76(3), 1373-1376.
- Williams, T. I., Combs, J. C., Thakur, A. P., Strobel, H. J., & Lynn, B. C. (2006). A novel Bicine running buffer system for doubled sodium dodecyl sulfate–polyacrylamide gel electrophoresis of membrane proteins. *Electrophoresis*, 27(14), 2984-2995.
- Wilson, W. W. (2003). Light scattering as a diagnostic for protein crystal growth—A practical approach. *Journal of Structural Biology*, 142(1), 56-65.
- Wiltfang, J., Arold, N., & Neuhoff, V. (1991). A new multiphasic buffer system for sodium dodecyl sulfate-polyacrylamide gel electrophoresis of proteins and peptides with molecular masses 100 000–1000, and their detection with picomolar sensitivity. *Electrophoresis*, 12(5), 352-366.
- Yamamoto, K., & Ogawa, K. (1983). Effects of NaOH-PIPES buffer used in aldehyde fixative on alkaline phosphatase activity in rat hepatocytes. *Histochemistry*, 77(3), 339-351.
- Yamanaka, T. (2003). The effect of pH on the growth of saprotrophic and ectomycorrhizal ammonia fungi in vitro. *Mycologia*, 95(4), 584-589.
- Yan, L. J., & Sohal, R. S. (1998). Mitochondrial adenine nucleotide translocase is modified oxidatively during aging. *Proceedings of the National Academy of Sciences*, 95(22), 12896-12901.
- Yang, X., & CHASTEEN, N. (1999). Ferroxidase activity of ferritin: effects of pH, buffer and Fe (II) and Fe (III) concentrations on Fe (II) autoxidation and ferroxidation. *Biochem. J.*, 338, 615-618.

- Yeoh, H. H., Badger, M. R., & Watson, L. (1981). Variations in kinetic properties of ribulose-1, 5-bisphosphate carboxylases among plants. *Plant Physiology*, 67(6), 1151-1155.
- Zaitseva, J., Holland, I. B., & Schmitt, L. (2004). The role of CAPS buffer in expanding the crystallization space of the nucleotide-binding domain of the ABC transporter haemolysin B from *Escherichia coli*. *Acta Crystallographica Section D: Biological Crystallography*, 60(6), 1076-1084.
- Zemann, A. J. (2001). Conductivity detection in capillary electrophoresis. *TrAC Trends in Analytical Chemistry*, 20(6), 346-354.
- Zhang, C., Baumgartner, R. A., Yamada, K., & Beaven, M. A. (1997). Mitogen-activated Protein (MAP) kinase regulates production of tumor necrosis factor- $\alpha$  and release of arachidonic acid in mast cells indications of communication between p38 and p42 Map kinases. *Journal of Biological Chemistry*, 272(20), 13397-13402.
- Zigler Jr, J. S., Lepe-Zuniga, J. L., Vistica, B., & Gery, I. (1985). Analysis of the cytotoxic effects of light-exposed HEPES-containing culture medium. *In Vitro Cellular & Developmental Biology*, 21(5), 282-287.